

# How to Install Freeze Dry Filter Paper

*During the freeze dry process, many samples may become a powder. Once this happens, the sample could be lost into the vacuum. In these instances, the filter paper that is shipped with the Fast-Freeze flasks can be used. The following steps show the correct way to insert the filter paper into the freeze dry flask lid.*

## STEP 1

*Place the lid on a flat surface.*



## STEP 2

*Place the filter paper over the opening of the lid.*



## STEP 3

*Push the filter paper into the lid using an adapter.*



## STEP 4

*Make sure the filter paper completely covers the hole once installed.*



## WARNING!

*DO NOT INSTALL  
THE FILTER PAPER  
ON THE UNDERSIDE  
OF THE LID.*



*Protecting your  
laboratory environment*

**LABCONCO**

**Labconco Corporation**

8811 Prospect Avenue • Kansas City, MO 64132-2696

816-333-8811 or 800-5525 • FAX: 816-363-0130

[www.labconco.com](http://www.labconco.com)