

LABCONCO CORPORATION 8811 Prospect Ave, Kansas City, MO 64132  
(816) 333-8811, Fax (816) 363-0130, (800) 821-5525

## **INSTALLATION INSTRUCTIONS FOR KIT #3697500 IV BAR KIT**

1. Before beginning, identify the kit(s) that you have selected. For installation in 3 and 4 foot wide Clean Benches (model numbers 36000xx or 36125xx. For 6 foot wide units (model number 36100xx), up to two IV bar kits can be installed.
2. Turn off the Clean Bench.
3. Install the IV Bar(s) by holding it in place in the rear of the work area near the diffuser. Use the four screws supplied in this kit to fasten the bar, using the predrilled holes in the underside of the canopy, as shown in Figure 1.

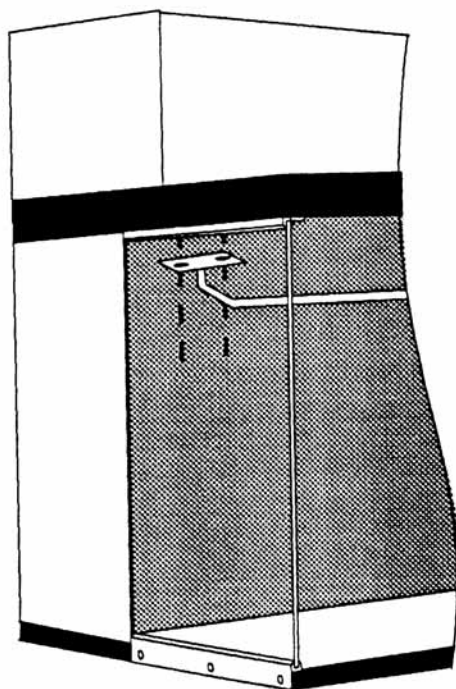


Figure 1