

Class II, Type C1 Biosafety Cabinet: Versatility, Safety & Savings Comparison

Abstract

Comparisons of Class II, Type A, Type B and Type C biosafety cabinets discern gaps in laboratory cost, safety, energy conservation and versatility. These four facets have become parameters by which all laboratory equipment is weighed and measured. The following gap analysis highlights the strengths and limitations of traditional Class II Type A and B biosafety cabinets (BSC's) and differentiates the role of the Type C1 BSC. The Type C1 bridges the gap between the recirculating Type A's and the vented Type B's and brings along new safety features and energy savings.



Introduction

Just when you thought you knew all aspects of the various types of Class II biosafety cabinets, along comes something truly innovative. Biosafety cabinets have experienced some general upgrades over recent years. Controls have become more interactive and energy-saving blowers have given us a reason to replace older cabinets. However, the core design of cabinet airflow has remained relatively unchanged.

However, Labconco introduced a new cabinet in 2014 that eases the selection process for Class II BSC's. The novel design drove industry experts to conclude that there are still new and enhanced ways to move air. Many industry experts agreed that this new cabinet fit neither the definition of a Type A, nor the definition of a Type B. Its unique airflow and design required it to have an entirely new classification – welcome the Class II, Type C1 biosafety cabinet (BSC).

This new Class II subtype's unique airflow configuration, installation, and exclusive ability to maintain safety even during emergency situations, compelled NSF International® to adopt, define and designate the new type of BSC in NSF/ANSI Standard 49. The product development team at Labconco Corporation describe the benefits of the Type C1 as, “unparalleled with conventional cabinets when summing up versatility, safety and savings. If you took the best features from A2s, B1s, B2s and put them all in a bowl with some new ideas and mixed them up you have the Type C1.”

New Concept in BSC Airflow

New technologies or re-adaptations of existing technologies break paradigms. Embracing the concept of the Type C1 BSC requires a paradigm shift.

From a distance, the Axiom® Class II, Type C1 takes on the appearance of other contemporary Labconco BSC's. What sets this cabinet apart is a patented “zoned” work area. The center two thirds of the work surface is clearly bordered by a pattern of intake slots. Called the Chem- Zone™, the center section is the dedicated exhaust portion of the cabinet

(Figure 1). Particulates and vapors produced in or near the Chem-Zone are ensured a direct path to the exhaust filter and out of the cabinet. Within this zone, there is no recirculation of air. Air outside this zone – downflow air at each end of the work surface and air taken in at the front grille – is recirculated to save energy in ducted applications.

This approach to cabinet airflow allows the center workspace to function as a typical B2 cabinet and provides an outer area that functions as an A2 cabinet. This provides a staging area for procedures as well as a protected place to keep equipment and supplies. One simply performs any tasks that generate volatiles near the center two-thirds of the cabinet. In installations where the Type C1 is ducted to the outside, the Chem-Zone has the identical performance one would expect from a Type B2 cabinet, but without the high air flow demands.

New Concept in BSC Airflow
“If you took the best features
from A2s, B1s, B2s and put
them all in a bowl with some
new ideas and mixed them up,
you have the Type C1.”

increases. A large remote blower is required because, over time, the pressure caused by a loading HEPA filter may increase to 4.0 inches water gauge or greater. In contrast, the Type C1 does not require a large dedicated remote blower, yet allows for the safe removal of chemicals in a single pass. How? Because the total exhaust area of the BSC is limited to the center portion of the work surface, exhaust air volume is reduced to approximately 40% of a comparable Type B2 BSC. Furthermore, a second blower, positioned downstream of the exhaust HEPA filter, pulls air through the BSC's exhaust HEPA filter, and delivers the air to the ductwork, where house exhaust can take over.

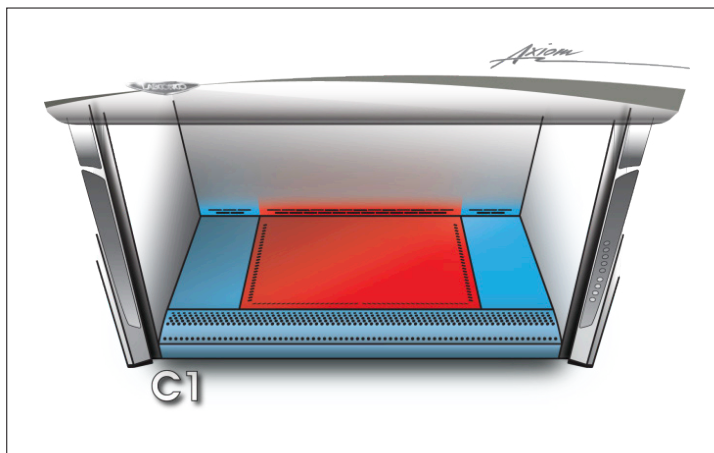


Figure 1 – The Chem-Zone™ (red) is a dedicated exhaust area of the work zone where vapor-forming chemicals are to be handled. The wings and grille (blue) provide supply air for the cabinet's sterile down flow supply air.

The Axiom® Type C1 challenges the thought that applications typically suitable for a Type B must have a “hard ducted,” dedicated exhaust run and dedicated exhaust fan. Traditional Type B biosafety cabinets require a hard duct (no canopy) because of the high vacuums necessary for their operation. On Type B BSC's, a dedicated roof-mounted blower, above the BSC, pulls the air through the BSC's exhaust HEPA filter. As the HEPA filter loads, the required vacuum

The Type C1 has caused institutions to rethink their requirements for exhausted biosafety cabinets. It has encouraged facility and lab managers reconsider which cabinet type they will want to own, whether it vents to the outside or not (Table 1).

Table 1 – Class II Gap Analysis.

Class II Type	A2	A2 with Canopy	B1	B2	A-mode	C1 B-mode
Exhausts to	Room	Outside	Outside	Outside	Room	Outside
Approximate exhausted volume (CFM)	N/A	270-400 CFM	270-400 CFM	660-800 CFM	N/A	300-400 CFM
Approximate vacuum req's (static pressure - in. w.g.)	N/A	0.15-0.25	0.7	1.6-4.0	N/A	0.3
Requires dedicated remote blower exhaust system	N/A	No	Yes	Yes	N/A	No
Volatile chemicals and radionuclides	No	Yes	Yes	Yes	No	Yes
Safe dedicated exhaust for chemicals	N/A	No	Yes**	Yes	N/A	Yes

This analysis looks at the impact each Class II biosafety cabinet subtype has on both safety and ventilation of the laboratory. Historically, a series of tradeoffs would have to be considered between safety, efficiency and expense. The new Type C1 dispels that notion, allowing a simple choice that offers safety, efficiency and cost savings, while adding flexibility.

*Data based on average of several 4' models for various subtypes.

** Type B1 BSCs do not have a clearly delineated or labeled dedicated exhaust area.

Versatility

Before the Type C1, deciding between a Type A and Type B required considerable forethought. Does one select the more economical Type A BSC or opt for the more costly Type B that offers a broader application potential? Making the right decision required answering questions such as: Will my researchers be using toxic volatile chemicals? Will I need to exhaust air out of the building for the life of this Type B BSC? If I commit to exhausting chemicals, will I require the installation of an additional exhaust duct and blower system? What will be the needs when this researcher moves out or changes applications? Since the Type C1 can operate as either a recirculating or exhausted cabinet, the buying decision is simplified and the risk of choosing the wrong cabinet eliminated. Here's why, the Type C1:

- Can easily convert from Type A mode (exhausting back into the laboratory) to Type B mode (venting exhaust to the outside), and vice versa.
- Can be connected to general laboratory exhaust systems, including those with fume hoods. Type B1 and B2 installations require dedicated remote exhaust systems for each cabinet.
- Can be connected to laboratory exhaust systems, the Type C1 can share ductwork with other equipment such as other BSC's or fume hoods.
- Can operate at the 8- or 10-inch sash opening. No other BSC can change sash heights in the field.

How does the Axiom achieve this versatility?

- Converting the Axiom from Type A mode to Type B mode is as “easy” as connecting or disconnecting the exhaust duct, selecting the exhaust preference in the software and having a certified technician verify inflow. The Type C1 BSC has the flexibility to change as the work in the lab changes.
- The reason for dedicated remote exhaust systems on Type B BSC's are their inherently large negative static pressures ($>1.5''$ W.C. or 0.37 kPa) along with airflow variations caused by other fume hoods or BSC's on the exhaust system. In contrast, the new Type C1 claims extremely low exhaust static pressures in the range of $0.2 - 0.4''$ W.C. (or < 0.1 kPa) while maintaining constant volume independent of the building exhaust.
- Unlike Type B1 and B2 cabinets, where personnel protective inflow velocities are directly influenced by fluctuations in the building exhaust, the Type C1 delivers a constant inflow regardless of drifts in the building's system. Additionally, it accomplishes this difficult task without the need for airflow or pressure sensors. Such devices require periodic calibration and replacement in order to keep the BSC operating as designed. (Hunter, et al., 2010).
- The Axiom, Class II, Type C1 has a software program selection that allows the transition from 8-inch to 10-inch operation to occur during any certification.

In review, these benefits allow the NSF listed Type C1 to function in Type B mode for maximum safety with volatiles or fumes and do so at a 10-inch sash height. Due to the exhaust requirements stated above, Type B cabinets only exist in 8-inch operational sash heights.

Unconventional Wisdom

First consider modern convention for Type B BSC's: personnel and product protection is entirely dependent upon the remote blower's proper operation. The ability for the Type B to protect the worker is challenged when there is a problem with the building exhaust system. There are a series of events that must happen quickly in the BSC for a positive user outcome. Current standards require ducted Type B cabinets to recognize when there is a loss of vacuum/flow (usually at less than 80% of exhaust capacity) and signal the cabinet to go into exhaust alarm within 15 seconds (NSF International, 2020). The design and calibration of the Type C1 allow that failure to be detected much sooner.

Centers for Disease Control and Prevention write that with Type B cabinets (U.S. Department of Health and Human Services, CDC, 2020):

For a Type B BSC, a significant loss of exhaust airflow will result in an alarm, turning off the cabinet blower(s). This stops the flow of air into the front of the BSC (inflow), negating personnel protection, potentially allowing materials in the total work area of the BSC to escape into the laboratory.

The momentum of the supply blower and the mass of downflow air may create a condition referred to as “puff back,” which is visualized during airflow visualization studies using smoke.

Labconco's solution to “puff-back” opposes conventional thinking of shutting down the cabinet as soon as possible during a building exhaust failure. Unlike a Type B2, the Type C1, equipped with its own exhaust blower, provides the means to maintain air flow; hence maintaining containment. Facility managers and safety officers must understand how the BSC reacts to loss of duct vacuum, and program the alarm conditions accordingly.

Solving “Puff-Back”

Puff-back occurs when a Class II, Type B cabinet’s remote exhaust system fails. As the internal blower windmills down, it continues to push air through the cabinet, and with no exhaust to remove that air, can result in a flow of air from the work area back into the laboratory. Even with the best BSC, the risk of “*puff-back*” can occur. To remedy this, Labconco developed a programmable active protection sequence. This system utilizes Constant Airflow Profile™ (CAP) to maintain safe inflow, downflow and exhaust through the cabinet. The safety officer is able to program how long this active protection is engaged between 0-300 seconds.

- IF THE TYPE C1 IS ALONE and not ganged (or manifolded) with other exhaust systems, then the full 5 minutes can be allotted to allow BSC operators ample time to cease work, perform emergency containment protocols and remove themselves from the cabinet safely.
- IF THE TYPE C1 IS GANGED with other ventilation or exhausted safety enclosures, then the appropriate period of time should be programmed allowing for the greatest protection duration at the face of the BSC while not violating safety protocols for other equipment.

A risk assessment is recommended to evaluate the possible outcomes in the event of building exhaust blower failure.

This argues against another commonly held rule in lab design and operations: a lab should never operate with positively pressurized fume hood or biosafety cabinet exhaust systems. Positive pressure can lead to chemicals being introduced to the building interior through openings or leaks. Following the loss of flow from a remote blower, Labconco realizes that the Type C1’s internal exhaust blower will, to some extent, positively pressurize the building duct for a short period of time. If the cabinet is connected to a shared duct system with other ventilation equipment, positive pressure can cause back-flows through other connected hoods; however, the pressure of the lab or room itself is more critical to determining the flow of air in these devices (U.S. Department of Health and Human Services, CDC, 2020; and Hunter, et al., 2016).

The recommended risk assessment will highlight the greatest potential for immediate harm. Is the most concentrated and immediate danger to life or health the loss of hazardous materials through the front of the biosafety cabinet? If so, then those users must be provided the necessary time to close containers and wipe surfaces. Operations being carried out in fume hoods during a building exhaust blower failure must be shut

The design and calibration of the Type C allow [a remote exhaust] failure to be detected much sooner.

down regardless of adjacent biosafety cabinets. The Type C1’s delayed shut down affords personnel time to manage the emergency situation systematically, control the most hazardous chemicals and biohazards first, and then carry out other shutdown protocols wearing the proper PPE.

Safety

The Labconco Axiom deliberately maintains the cabinet's exhaust blower for 10 seconds after the supply downflow blower has shut down. This patent-pending feature ensures the interior remains negatively pressurized and eliminates any possibility for "puff-back." Additionally, the Type C1 has a programmable shutdown timer. The countdown timer should be set to the shortest possible interval required for lab personnel to secure hazards inside the cabinet upon an emergency.

To be clear, the Type C1 is not designed to operate indefinitely with a failed building exhaust blower. The BSC will not start up until it detects adequate exhaust; furthermore, the maximum period of time that the cabinet can be programmed to maintain containment during an exhaust alarm is five minutes. When the shutdown timer reaches zero, the cabinet will sequentially turn off the supply blower followed by the cabinet's exhaust blower 10 seconds later. The cabinet will not return to operation until it is restarted by the operator and the building exhaust system problem has been remedied.

All biosafety cabinets that have met the standards of NSF/ANSI 49 or EN 12469 and have been properly certified by a qualified technician should safely protect the operator and the work samples. The question is where can new biosafety cabinet technology exceed the accepted norm? Labconco's Type C1 advances BSC safety in two ways. First, by enhancing the handling of emergency situations and secondly by ensuring consistent airflow performance between periodic validations by certifiers.

- The Labconco Axiom BSC is able to maintain complete containment and product protection upon exhaust system failure. Instead of attempting to shut down the cabinet quickly as all other Type B cabinets do, it does the opposite.
- More advanced controls available in the Axiom essentially eliminates adjustments once required of certifiers.

The Axiom makes necessary adjustments to inflow and downflow to maintain normal operating airflows. The lab personnel are instantly made aware of the building exhaust failure audibly and visually through the line-of-sight color display. A highly visible countdown timer affords the researcher a preset 0 to 300 seconds to secure hazardous or valuable materials and surface decontaminate before the cabinet shuts down.

Furthermore, the Axiom carries on the proven reliability of earlier Labconco biosafety cabinets with high quality ECM (electronically commutated motors) driven blowers. The same Constant Airflow Profile™ (or CAP) first utilized in 2007 on the Logic® (Hunter, et al., 2008) ensures the blowers deliver the same volume of air despite blockages in the grille or HEPA/ULPA filter loading. The inflow and downflow volumes remain constant between periodic certification inspections.

Savings

How would the Type C1 compare in total cost of ownership with conventional biosafety cabinet types? Since it has the ability to exhaust back into the laboratory like an A2, or the adaptability to be ducted for single-pass exhaust from the center of the work surface like a B2, let's view it both ways.

Certainly in terms of electrical power usage and the loss of conditioned air, the most energy efficient way to operate a BSC is to return the exhaust air back into the laboratory. Returning the air to the room is not acceptable when the application involves volatile toxic chemicals, odors or radionuclides.

When the cabinet is exhausting back into the room, the Type C1 consumes the same energy as any of the energy-efficient A2 cabinets that utilize ECM blower motors.

Energy savings is readily apparent in the Type C1 when used in an application that requires remote exhaust. In terms of removing conditioned air from the building, the Type C1 exhausts similar air volumes as do Type B1s and A2s. These cabinets remove 50% less air than does a Type B2. Employing any one of them will reduce operating cost by half. (Table 2)

The “green” design intent for the Type C1 is to permit the facility manager along with the research managers to optimize energy savings and allow the Type C1 to run in ‘Type A mode’. This is done by disconnecting Type C1 cabinets from exhaust ducts where possible. Applications change; why not employ a biosafety cabinet that can adapt?

Further comparison with Type B1 or Type B2 cabinets reveals the potential for greatly reduced expenses in design and construction. The Type C1 is capable of being connected to typical ganged laboratory exhaust systems thereby eliminating the need for additional dedicated duct and blower systems. The savings comprised of consulting services, utility space, heavy ducting, exhaust blower, installation, wiring and roof penetrations are substantial.

Table 2 – Life Cycle Cost Analysis.

Class II Type	A2	A2 with Canopy	B1	B2	C1	
					A-mode	B-mode
Upfront cost of installation*	\$300	\$400	\$5,150	\$5,150	\$300	\$400
Lifetime maintenance Cost (Service & HEPA/ULPA filters)	\$4,500	\$4,500	\$4,500	\$4,500	\$4,500	\$4,500
Lifetime operating cost (based on \$8 / CFM / year)	N/A	\$40,500	\$40,500	\$87,000	N/A	\$42,000
ESTIMATED TOTAL COST (INSTALLATION & OPERATION)	\$4,800	\$45,400	\$50,150	\$96,650	\$4,800	\$46,900

With purchase price and cost an ever mounting concern for capital laboratory equipment, it is important to know what the different Class II subtypes will cost through the life of the product (based on the 15 year life of a 4' BSC, 20' of epoxy-coated duct for exhausting out of a single story). Costs are approximate only.

*Cost of installation includes materials for ductwork, exhaust stack, and labor on a single cabinet.

Conclusion

Occasionally products emerge that have the potential to provide an option not previously available in our industry. We believe the Labconco Axiom, Class II, Type C1 is the new generation of BSC in a world that is ready to evolve. New versatility allows the BSC to replicate the strengths of the Type A and Type B cabinets all while addressing each of their limitations. The Axiom also provides lifelong flexibility to convert at any time

necessary. The cost of ownership is far less than conventional Type B BSC's. Low exhaust volume, the lowest required vacuums and the ability to connect to ganged exhaust systems add up to big savings.

Finally, the purpose for biosafety cabinets is safety. Original thinking and new technology provide consistent containment and product protection even in emergency situations. The Type C1 is a genuine advancement in bio-containment that deserves attention.

Works Cited

Garrett, B., Labconco Corporation. (2013). *Maximizing User Safety Through Human Factors Design* [White paper]. Retrieved from <https://www.labconco.com/resources/white-papers>.

Hunter, J., Hays, B. & Rouse, M. (2016). *Effect of the Type C1 Active Protection Protocol on a Disabled Exhaust System and Other Proximal Ventilation Devices*. Kansas City, MO: Labconco Corporation.

Hunter, J., Meinders, M. & Garrett, B. (2010). *Controlling Airflow in Class II Biosafety Cabinets*. Kansas City, MO: Labconco Corporation.

Hunter, J. & Rouse, M. (2008). *The Advantages of Using an Electronically Commutated Motor (ECM) in Biosafety Cabinets*. Kansas City, MO: Labconco Corporation.

NSF International. (2021). *Biosafety Cabinetry: Design, Construction, Performance and Field Certification* (NSF/ANSI 49 - 2020). NSF International.

U.S. Department of Health and Human Services/Center for Disease Control. (2020). *Biosafety in Microbiological and Biomedical Laboratories*. 6th ed. Washington, D.C.: U.S. Government Printing Office.

LABCONCO CORPORATION

8811 Prospect Avenue
Kansas City, MO 64132
(800) 821-5525 | (816) 333-8811
labconco.com



©2022 LABCONCO CORPORATION
Printed in the U.S.A.
Design subject to change without notice.
20221110