LABCONCO CORPORATION

8811 Prospect Avenue
Kansas City, MO 64132
(800) 821-5525, (816) 333-8811
(816) 363-0130 fax
labconco@labconco.com

User's Manual

Protector[®] PVC Laboratory Fume Hoods

Perchloric Acid Models

14041 Series
14051 Series
14061 Series
14081 Series

Acid Digestion Models

14141 Series
14151 Series
14161 Series
14181 Series

To receive important product updates,
complete your product registration card
online at register.labconco.com

Please read the User's Manual before operating the equipment.

Copyright © 2021 Labconco Corporation. The information contained in this manual and the accompanying products are copyrighted and all rights reserved by Labconco Corporation. Labconco Corporation reserves the right to make periodic design changes without obligation to notify any person or entity of such change.

Warranty

Labconco Corporation provides a warranty to the original buyer for the repair or replacement of parts and reasonable labor as a result of normal and proper use of the equipment with compatible chemicals. Broken glassware and maintenance items, such as filters, gaskets, light bulbs, finishes and lubrication are not warranted. Excluded from warranty are products with improper installation, erratic electrical or utility supply, unauthorized repair and products used with incompatible chemicals.

The warranty for Protector® PVC Laboratory Fume Hoods will expire one year from date of installation or two years from date of shipment from Labconco, whichever is sooner. Warranty is non-transferable and only applies to the owner (organization) of record.

Buyer is exclusively responsible for the set-up, installation, verification, decontamination or calibration of equipment. This limited warranty covers parts and labor, but not transportation and insurance charges. If the failure is determined to be covered under this warranty, the dealer or Labconco Corporation will authorize repair or replacement of all defective parts to restore the unit to operation. Repairs may be completed by 3rd party service agents approved by Labconco Corporation. Labconco Corporation reserves the rights to limit this warranty based on a service agent's travel, working hours, the site's entry restrictions and unobstructed access to serviceable components of the product.

Under no circumstances shall Labconco Corporation be liable for indirect, consequential, or special damages of any kind. This warranty is exclusive and in lieu of all other warranties whether oral, or implied.

Returned or Damaged Goods

Do not return goods without the prior authorization from Labconco. Unauthorized returns will not be accepted. If your shipment was damaged in transit, you must file a claim directly with the freight carrier. Labconco Corporation and its dealers are not responsible for shipping damages.

The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

Limitation of Liability

The disposal and/or emission of substances used in connection with this equipment may be governed by various federal, state, or local regulations. All users of this equipment are required to become familiar with any regulations that apply in the user's area concerning the dumping of waste materials in or upon water, land, or air and to comply with such regulations. Labconco Corporation is held harmless with respect to user's compliance with such regulations.

Contacting Labconco Corporation

If you have questions that are not addressed in this manual, or if you need technical assistance, contact Labconco's Customer Service Department or Labconco's Product Service Department at 1-800-821-5525 or 1-816-333-8811, between the hours of 7:30 a.m. and 5:30 p.m., Central Standard Time.

TABLE OF CONTENTS

CHAPTER 1: INTRODUCTION	1
About This Manual	2
Typographical Conventions	3
(Conventions Typographiques)	
Your Next Step	5
CHAPTER 2: PREREQUISITES	6
Location Requirements	7
Support Requirements	7
Exhaust Requirements	7
Exhaust Washdown Requirements	8
Electrical Requirements	9
(Exigences électriques)	
Service Line Requirements	9
Space Requirements	10
CHAPTER 3: GETTING STARTED	11
Unpacking Your Fume Hood	12
Removing the Shipping Skid	12
Sash Weight Release	13
Installing the Hood on a Supporting Structure & Hardboard Work Top	13
Connecting to the Hood Exhaust System	16
Connecting to the Exhaust Washdown	17
Connection to the Washdown Drain	17
Connecting the Electrical Supply Source	18
Connecting the Service Lines to the Protector Hood	20
Sealing the Protector Hood to the Work Surface	21
Certifying the Protector Fume Hood	21
CHAPTER 4: PERFORMANCE FEATURES AND SAFETY	
PRECAUTIONS	22
Performance Features	22
Safety Precautions	27
(Consignes de sécurité)	

CHAPTER 5: USING THE PROTECTOR PVC	32
Operating the Vertical-Rising Glass Sash	32
Operating the Vertical-Rising Lexan Sash	32
Operating the Blower	32
Operating the Lights	32
Operating the Washdown Control Valve	33
Working in Your Protector PVC Fume Hood	33
CHAPTER 6: MAINTAINING THE PROTECTOR PVC	35
Service Safety Precautions	35
(Mesures de sécurité des services)	
Routine Maintenance Schedule	36
Routine Service Operations	37
CHAPTER 7: MODIFYING THE PROTECTOR PVC	38
Installing Hardboard Support Surface	38
Installing PVC Eco Foil Made from PVC Material	38
Installing Ceiling Enclosures Above the Fume Hood	39
Installing Rear Panels Behind the Fume Hood	41
Installing Guardian Airflow Monitors	41
Sash Stop Kit – Field Installation	42
Installing an Electrical Duplex Outlet	42
(Installation d'une sortie Duplex électrique)	
Installing the Acid Vent Kit from the Acid Cabinet	42
Installing Additional Washdown Valve	42
CHAPTER 8: TROUBLESHOOTING	43
APPENDIX A: PROTECTOR PVC COMPONENTS	46
APPENDIX B: PROTECTOR PVC DIMENSIONS	49
APPENDIX C: PROTECTOR PVC SPECIFICATIONS	53
(ANNEXE C: SPÉCIFICATIONS PROTECTEUR DU PVC MINISTRE)	
APPENDIX D: SERIAL NUMBER TAG DESCRIPTION	56
(ANNEXE D: DESCRIPTION DE SÉRIE TAG NUMBER)	
APPENDIX E: REFERENCES	62
APPENDIX F: VAV OPEN AREA GUIDELINES	65

CHAPTER 1

INTRODUCTION

Congratulations on your purchase of a Labconco Protector® PVC Perchloric Acid Laboratory Fume Hood or Protector® PVC Acid Digestion Laboratory Hood. The difference between the two hoods is that the Protector PVC Perchloric Acid Hood uses a glass sash and the Protector PVC Acid Digestion Hood uses a Lexan™ sash to combat the fogging of glass caused by hydrofluoric acid (HF). Note that Perchloric acid will react with the Lexan sash and is not to be used in an Acid Digestion Hood. The Protector PVC Perchloric Laboratory Hood is designed to efficiently ventilate procedures using Perchloric acid. The Protector PVC Acid Digestion Laboratory Hood is designed to efficiently ventilate all other heavy acid operations including hydrofluoric acid. It is the result of Labconco's more than 50 years experience in manufacturing fume hoods, and users like you suggested many of its features to us.

The Protector PVC Perchloric Acid Fume Hood has been engineered to provide maximum safety in a laboratory while using Perchloric acid. **Consult your safety officer for appropriate chemical combinations with Perchloric acid. Perchloric acid is a strong oxidizing agent, and should not be used in combination with any organic material as the possibility of a hazardous chemical reaction or explosion could develop.** The Protector PVC Acid Digestion Fume Hood has been designed to handle most other types of acids including hydrofluoric acid (HF), but not Perchloric acid. The Protector PVC Hoods offer many unique features to enhance safety, performance, and visibility.

To take full advantage of them, please acquaint yourself with this manual and keep it handy for future reference. If you are unfamiliar with how fume hoods operate, please review *Chapter 4: Performance Features and Safety Precautions* before you begin working in the fume hood. Even if you are an experienced fume hood user, please review *Chapter 5: Using The Protector PVC*, which describes the Protector PVC Perchloric Acid Hood features so that you can use the hood efficiently.

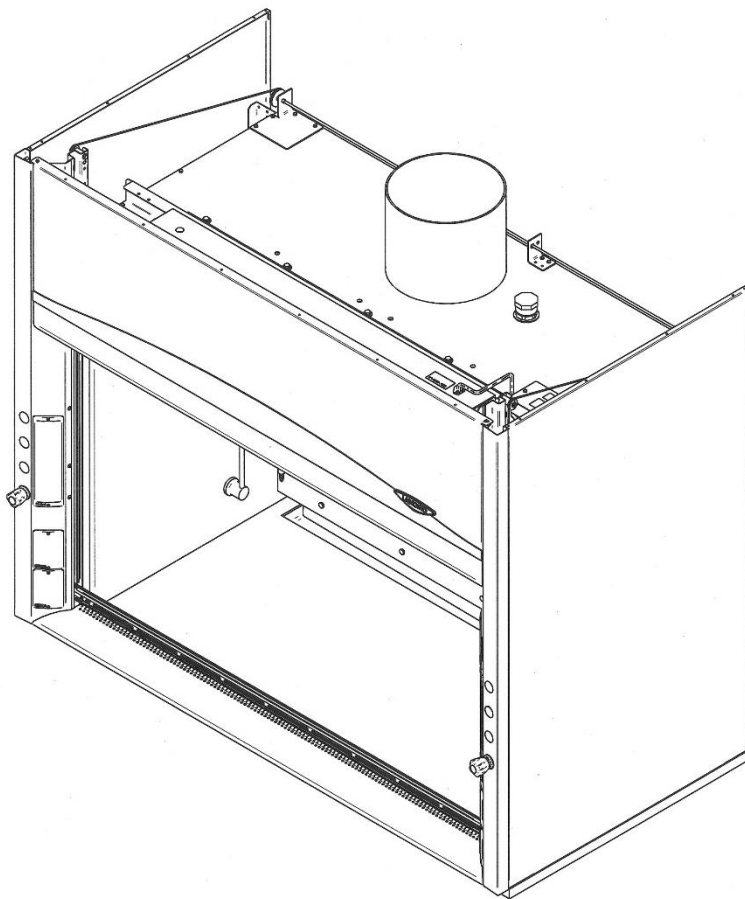
Figure 1-1

About This Manual

This manual is designed to help you learn how to install, use, and maintain your laboratory fume hood. Instructions for installing optional equipment on your hood are also included.

Chapter 1: Introduction provides a brief overview of the laboratory fume hood, explains the organization of the manual, and defines the typographical conventions used in the manual.

Chapter 2: Prerequisites explains what you need to do to prepare your site before you install your laboratory fume hood. Electrical and service requirements are discussed.

Chapter 3: Getting Started contains the information you need to properly unpack, inspect, install, and certify your laboratory fume hood.

Chapter 4: Performance Features and Safety Precautions explains how the Protector operates and the appropriate precautions you should take when using the fume hood.

Chapter 5: Using The Protector PVC discusses the basic operation of your fume hood. Information on how to prepare, use and shut down your Protector Hood are included.

Chapter 6: Maintaining The Protector PVC explains how to perform routine maintenance on your fume hood.

Chapter 7: Modifying The Protector PVC explains how to modify the fume hood or add accessories.

Chapter 8: Troubleshooting contains a table of problems you may encounter while using your laboratory fume hood including the probable causes of the problems and suggested corrective actions.

Appendix A: Protector PVC Components contains labeled diagrams of all of the components of the fume hoods.

Appendix B: Protector PVC Dimensions contains comprehensive diagrams showing all of the dimensions for the laboratory fume hoods.

Appendix C: Protector PVC Specifications contains the electrical requirements for laboratory fume hood. Wiring diagrams are also included.

Appendix D: Serial Number Tag Description provides current rating code used on serial number tag.

Appendix E: References lists the various resources available that deal with laboratory fume hoods.

Appendix F: VAV Open Area Guidelines to properly assess fume hood effective area whenever variable air volume (VAV) mechanical systems are installed.

Typographical Conventions

Recognizing the following typographical conventions will help you understand and use this manual:

- Book, chapter, and section titles are shown in italic type (e.g., *Chapter 3: Getting Started*).
- Steps required to perform a task are presented in a numbered format.
- Comments located in the margins provide suggestions, reminders, and references.

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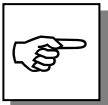
- Important information is presented in capitalized type in paragraphs that are preceded by the pointer icon. It is imperative that the information contained in these paragraphs be thoroughly read and understood by the user.

- The S icon indicates the text is specific to the standard model.

- The A icon indicates the text is specific to the A-Style Combination Sash Model.

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- CAUTION – Hot Surface.
- AVERTIR – Surface Chaude

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Your Next Step

If your Fume Hood needs to be installed, proceed to *Chapter 2: Prerequisites* to ensure your installation site meets all of the requirements. Then, go to *Chapter 3: Getting Started* for instructions on how to install your laboratory fume hood and make all of the necessary connections.

If you would like to review how laboratory fume hoods operate, go to *Chapter 4: Performance Features and Safety Precautions*.

For information on the operational characteristics of your laboratory fume hood, go to *Chapter 5: Using The Protector PVC*.

If your laboratory fume hood is installed and you need to perform routine maintenance on the cabinet, proceed to *Chapter 6: Maintaining Your Protector PVC*.

For information on making modifications to the configuration of your fume hood, go to *Chapter 7: Modifying The Protector PVC*.

Refer to *Chapter 8: Troubleshooting* if you are experiencing problems with your fume hood.

CHAPTER 2

PREREQUISITES

Before you install your laboratory fume hood, you need to prepare your site for installation. Carefully examine the location where you intend to install your hood. You must be certain that the area is level and of solid construction. In addition, a dedicated source of electrical power must be located near the installation site.

Carefully read this chapter to learn the requirements for your installation site:

- The location requirements.
- The support requirements.
- The exhaust requirements.
- The exhaust washdown requirements.
- The electrical power requirements.
- The service line requirements.
- The space requirements.

Refer to *Appendix B: Protector PVC Dimensions* for complete fume hood dimensions.

Refer to *Appendix C: Protector PVC Specifications* for complete laboratory fume hood electrical and environmental conditions, specifications and requirements.

Location Requirements

The fume hood should be located away from traffic patterns, doors, windows, fans, ventilation registers, and any other air-handling device that could disrupt its airflow patterns. All windows in the room should be closed.

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Support Requirements

DO NOT install the fume hood on a cart, dolly, or mobile bench. **ALL** Protector Hood installations must be permanent and stationary. The supporting structure usually consists of a base cabinet and chemically resistant work surface.

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Exhaust Requirements

The exhaust duct connection has been designed for 12" nominal duct (12.75" OD) to allow for minimum static pressure loss while operating at 100 fpm face velocities. The 12" diameter exhaust duct also allows for proper transport velocities away from the hood in the 1000 fpm to 2500 fpm range. The proper exhaust volume and static pressure loss are listed next for each hood model:

Chapter 2: Prerequisites

Airflows and Static Pressure

Hood Type	Face Velocity (fpm)		Eco Energy Saving Air Foil			
	Sash at	Sash at 62.5%	Airflow Volumetric Rate (CFM) @			
	Full open at 28"	open at 18"	Static Pressure (inches of water)			
			4' Hood	5' Hood	6' Hood	8' Hood
PVC Hoods						
Perchloric	125	200	905, 0.27"	1195, 0.44"	1475, 0.59"	2050, 0.35"
and	100	160	725, 0.17"	955, 0.28"	1180, 0.38"	1640, 0.22"
Acid	N/A	125	565, 0.11"	745, 0.17"	920, 0.23"	1280, 0.13"
Digestion	N/A	100	455, 0.07"	595, 0.11"	740, 0.15"	1025, 0.09"

12.75" OD

Exhaust Recommendations highlighted

Proper PVC blower selection can be determined from these exhaust requirements and the total system static pressure loss. Contact Labconco Customer Service for assistance in sizing a blower system. Refer to *Appendix F: VAV Open Area Guidelines* to properly assess the fume hood area opening, allowing the variable air volume (VAV) mechanical system to be programmed at various sash positions.

Exhaust Washdown Requirements

The entire exhaust duct system **must** feature an adequate washdown system for **Perchloric acid** applications, but are not always required (but recommended) for standard mineral acid applications such as hydrochloric, nitric or sulfuric. This includes the exhaust ductwork, blower, hood structure itself, and any ductwork located outside the blower. Wash rings are required to be placed in the exhaust duct system every 8 to 10 feet and after any change in direction to provide adequate washdown. Contact Labconco for washdown ring ordering information. Order part number 4746000 for a 10" wash ring and 4746100 for a 12" wash ring. Depending on the length of duct run and number of wash rings and elevation differences, Labconco recommends one additional washdown valve for every three wash rings. The washdown nozzles are designed at a normal pressure of 40 psi with a flowrate of 1 gallon per minute and a minimum pressure of 10 psi with a flow rate of 0.5 gallons per minute. The washdown nozzle at a maximum pressure of 100 psi has a flowrate of 1.5 gallons per minute. Each washdown valve can only produce up to 3.5 gallons per minute at 40 psi and additional washdown valves are ordered frequently and installed on the fume hood front corner post. Order 9808310 Washdown Valve Kit (3/8 lines) per sizing chart that follows.

Additional Washdown Valve Sizing Chart

Duct Run	<u>4' PVC</u>		<u>5' PVC</u>		<u>6' PVC</u>		<u>8' PVC</u>		
	<u>30'</u>	<u>60'</u>	<u>30'</u>	<u>60'</u>	<u>30'</u>	<u>60'</u>	<u>30'</u>	<u>60'</u>	
Nozzles Mounted on Fume Hood	3	3	4	4	4	4	6	6	
Additional Wash Rings	3	6	3	6	3	6	3	6	
Additional Washdown Valve Kits Required (9808310)							1	2	Minimum 10 psi pressure
	1	2	1	2	1	2	or	or	
							2	3	Normal 40 psi pressure

A washdown and drain should be provided in your exhaust blower. This washdown water and effluent material can be drained from the exhaust blower back into the exhaust duct for disposal. Contact Labconco for PVC blowers. The Protector PVC Hood washdown drain outlet is sized for 2" nominal (2.375" OD) PVC pipe.

Electrical Requirements

The Protector Hood models feature internal wiring for the LED light assembly and light switch. All internal wiring is terminated at the single point wiring junction box for hook-up by a qualified electrician. The blower switch, and light switch wires are also terminated at the single point wiring junction box for hook-up by a qualified electrician. Refer to *Chapter 3: Getting Started* and *Appendix C: Protector PVC Specifications* for the wiring diagram for proper electrical installation.

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Service Line Requirements

All service lines to the laboratory fume hood should be ¼ inch outside diameter, copper (brass for natural gas), and equipped with an easily accessible shut-off valve, should disconnection be required. Recommended operating pressure is 40 PSI, with a maximum allowable pressure of 200 PSI. Consider a pressure regulator to reduce line pressure to 40 PSI. Please check with local codes for other requirements.

Space Requirements

The dimensions for the different models are shown in *Appendix B: Protector PVC Dimensions*.

CHAPTER 3

GETTING STARTED

Now that the site for your laboratory fume hood is properly prepared, you are ready to unpack, inspect, install, and certify your unit. Read this chapter to learn how to:

- Unpack and move your Protector PVC Hood.
- Set up the fume hood with the supporting structure and hardboard work top.
- Connect to an exhaust system suitable for Perchloric Acid and acids.
- Connect the exhaust washdown to the exhaust system.
- Connect to the washdown drain.
- Connect the electrical supply source.
- Connect the service lines.
- Arrange certification of your Protector PVC Hood.

Depending upon which model you are installing, you may need common plumbing and electrical installation tools in addition to 5/16", 3/8", 7/16", and 1/2" wrenches, ratchets, sockets, a nut driver set, a flat-blade screwdriver, a Phillips screwdriver, and a carpenter level to complete the instructions in the chapter.

The Protector PVC Hood models weigh between 400 to 800 lbs. (182-363 kg). The shipping skid allows for lifting with a mechanical lift truck or floor jack. If you must lift the fume hood manually, follow safe-lifting guidelines. Normally, the fume hood can be slid off a hydraulic lift table and be placed into position on top of the work surface. Do not lift by the front air foil.

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

Unpacking Your Fume Hood

Carefully remove the shrink-wrap or carton on your fume hood and inspect it for damage that may have occurred in transit. If your unit is damaged, notify the delivery carrier immediately and retain the entire shipment intact for inspection by the carrier.

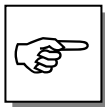The United States Interstate Commerce Commission rules require that claims be filed with the delivery carrier within fifteen (15) days of delivery.

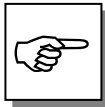DO NOT RETURN GOODS WITHOUT THE PRIOR AUTHORIZATION OF LABCONCO. UNAUTHORIZED RETURNS WILL NOT BE ACCEPTED.

IF YOUR HOOD WAS DAMAGED IN TRANSIT, YOU MUST FILE A CLAIM DIRECTLY WITH THE FREIGHT CARRIER. LABCONCO CORPORATION AND ITS DEALERS ARE NOT RESPONSIBLE FOR SHIPPING DAMAGES.

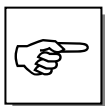
Do not discard the shipping skid or packing material for your fume hood until you have checked all of the components and installed and tested the unit. **The PVC fume hood baffle is installed as shown in Figure 1-1.** Do not remove the fume hood from its shipping skid until it is ready to be placed into its final location. Move the unit by placing a flat, low dolly under the shipping skid, or by using a floor jack.

Do not move the hood by tilting it onto a hand truck.

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Removing the Shipping Skid

LEAVE THE FUME HOOD ATTACHED TO ITS SHIPPING SKID UNTIL IT IS AS CLOSE TO ITS FINAL LOCATION AS POSSIBLE. MOVE THE HOOD BY USING A SUITABLE FLOOR JACK, OR BY PLACING A FURNITURE DOLLY UNDERNEATH THE SKID. DO NOT MOVE THE HOOD BY TILTING IT ONTO A HAND TRUCK.

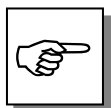
After you verify the fume hood components, move your hood to the location where you want to install it. Follow the steps listed next to remove the shipping skid from your unit.

1. Remove the side panels by unscrewing the Phillips screws.

2. Find the hardware (bolts, washers, nuts) that attach the fume hood to the skid and remove the hardware. Some hardware is on the sides and some is on the back.

Sash Weight Release

To protect the fume hood from damage in shipment, the sash weight has been secured to the back of the fume hood with screws. Simply remove the screws and make sure the sash cables or chains are on the pulleys or sprockets before operation of the sash.

NOTE: THE SASH WEIGHT ITSELF WAS INDIVIDUALLY MATCHED FOR THIS SPECIFIC HOOD AND SHOULD NOT BE EXCHANGED ON ANY OTHER UNIT.

Installing the Hood on a Supporting Structure and Hardboard Work Top

The Protector Hood is heavy! Use caution when lifting or moving the unit.

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When installing the Protector PVC Hood onto the lower hardboard work top, ensure that the structure can safely support the combined weight of the fume hood and any related equipment. **The hardboard work top is aligned flush with the back of the fume hood for good airflow: this will provide the correct spacing under the air foil for proper bypass airflow. The lower base cabinets are placed flush with the front of the hardboard work top as shown in Figure 3-2. The hardboard work top supports the PVC Hood and has a cutout to clear the drain trough.**

WARNING: It is important to support the rear of the work surface and fume hood. The cross support provides support for the bottom of the work surface. Install the cross support after the base cabinets and work surface are leveled and before installing the hood.

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Chapter 3: Getting Started

The following are instructions for mounting a cross support:

1. Level the base cabinets and the hardboard work top. Work top should be placed flush with the back of the fume hood as shown in Figure 3-1.
2. Scribe a line on the wall or back of the base cabinet to locate the support under the work surface.
3. Mount the support by attaching it to the wall or base cabinet.
4. Anchor cabinet(s) with seismic restraints if required by the local code in your area.

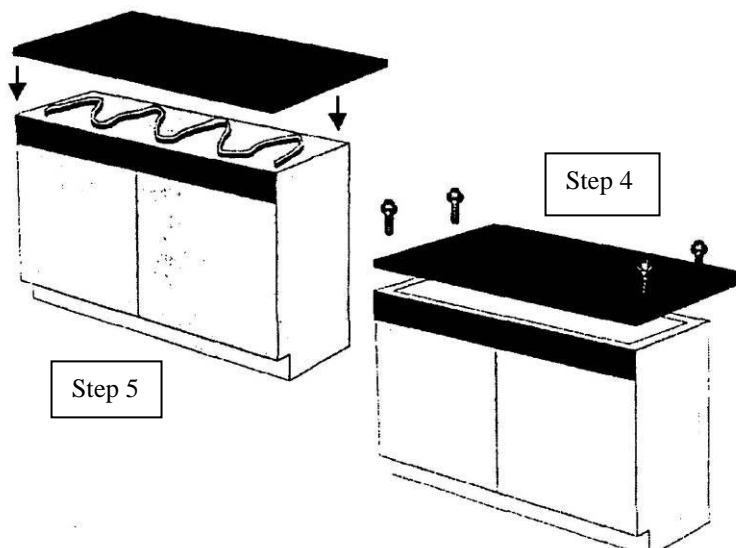
Figure 3-1

5. Secure the hardboard surface to the top of the base cabinet(s) using adhesives with the Acid Storage Cabinet(s). Be sure the PVC Hood drain trough is accounted for in the relief area of the hardboard top. See figure 3-2.
6. After completing step 5, apply a sufficient quantity of adhesive to the top surface of the hardboard.
7. Place the hood on top of the work surface and cross support.
8. Place heavy weights on the inside of the fume hood for approximately 24 hours. This provides a proper bond of the adhesive. Protect the hood's integral work surface from possible damage by wrapping any weighted material first.
9. Place the hood on top of the work top and cross support.

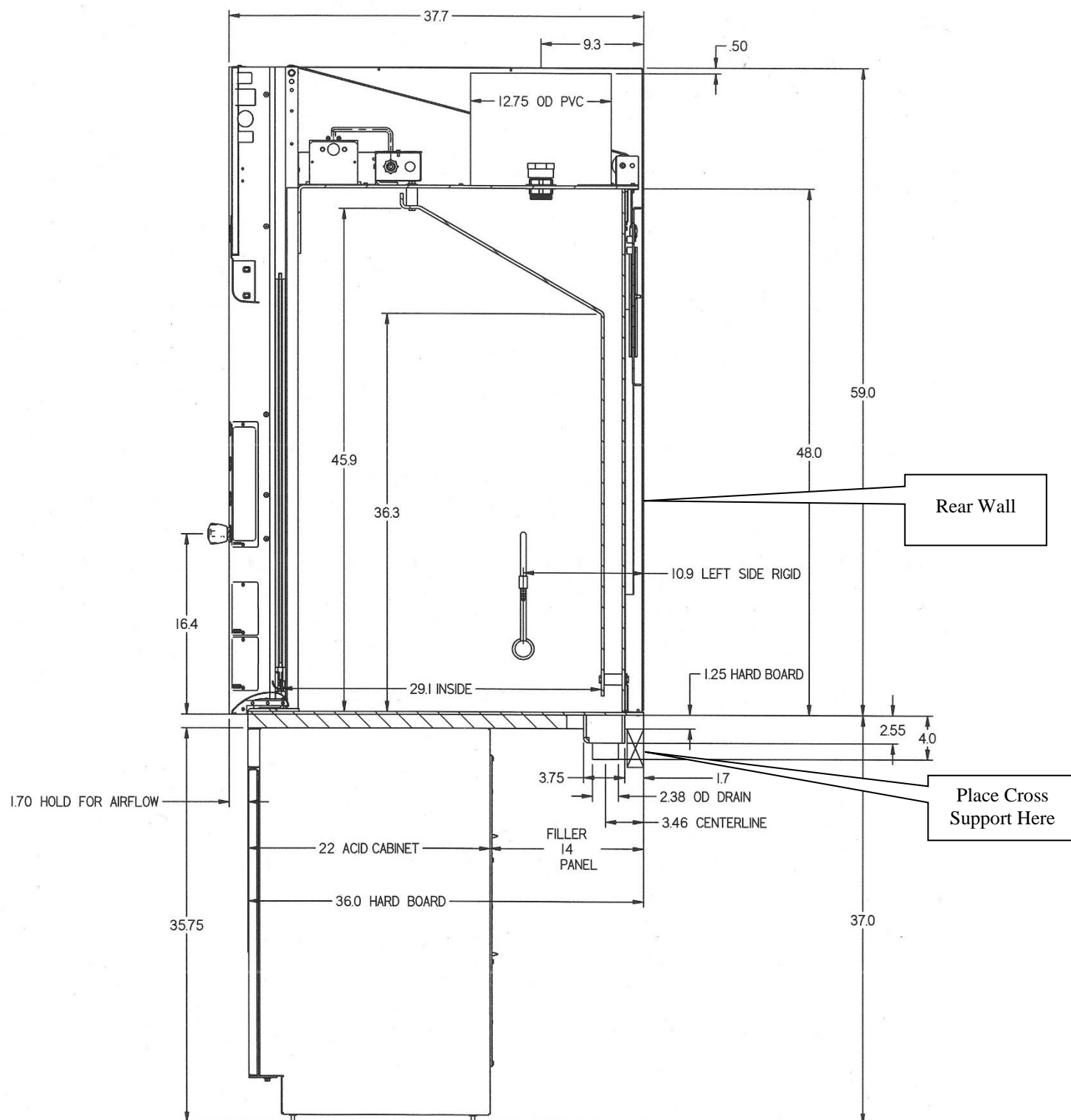
Figure 3-2

Connecting to the Hood Exhaust System

WARNING: The weight of the exhaust ductwork system must be supported independently of the hood superstructure. Do not allow this weight to be supported by the hood structure as damage to the hood may occur.

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The exhaust connection should be installed by a qualified HVAC contractor. The exhaust connection on your hood is manufactured with 12" PVC nominal pipe (12.75" OD) to allow for minimum static pressure loss with proper transport velocities away from the hood. Consult Labconco Customer Service should you require help sizing your blower for the exhaust volume and total system static pressure loss.

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The selected exhaust duct material must be fabricated out of type 1 unplasticized polyvinyl chloride or type 316 stainless steel material. The duct should be fully welded in the case of the stainless steel material or solvent bonded in the case of polyvinyl chloride so that it does not have any cracks or holes where Perchloric acid salts or acids can collect.

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The exhaust duct system for a PVC Fume Hood should be run in a vertical position and feature horizontal runs only where it cannot be avoided. Should a horizontal run be a part of your exhaust system, it should be sloped downward toward the hood to prevent the pooling of the washdown water and perchloric acid salts or acids in the duct.

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CAUTION: PVC Fume Hoods must never be manifolded into a common exhaust system. They require a dedicated exhaust system.

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Connecting to the Exhaust Washdown

The entire exhaust duct system must have an adequate washdown system. This includes the exhaust ductwork located between the exhaust blower and the hood structure, the exhaust blower, and any ductwork located on the outlet side of the blower. Wash rings are recommended to be placed in the exhaust duct system approximately every 8 to 10 feet and after every change in direction to provide an adequate washdown. A drain should be provided in your exhaust blower and this washdown water and effluent material can be drained from the exhaust blower back directly into the exhaust duct for disposal. The PVC Hood was designed with an auxiliary port plugged at the factory, and it is to be connected to the washdown valve system to provide water to the wash rings. The plug is located on top of the hood near the first washdown nozzle. Plumb from the plugged connector with a 3/8" rigid polyethylene, copper, or stainless steel line.

Connecting to the Washdown Drain

The Protector PVC Hood has been supplied with a 2" nominal (2.375" OD) PVC pipe drain. Connect the 2" pipe to your floor drain. Be sure the connection is either fully solvent bonded or leak tight. See Figure 3-2.

Connecting the Electrical Supply Source to the Protector Fume Hood

Prior to connecting any electrical wiring to the fume hood structure, refer to the hood identification plate for the proper electrical requirements of your specific model.

WARNING: The building electrical supply system for Protector Hoods is required to include overload protection. A switch or circuit breaker shall be in close proximity to the equipment and within easy reach of the operator. The switch or circuit breaker is to be marked as the disconnecting device for the equipment.

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The identification plate, model number, serial number, and electrical connection boxes are accessible from the front of the fume hood by removing the front panel.

The Protector PVC Hood is normally wired for 115 Volt, 50/60 Hz, 20 Amp or 230Volt, 50/60 Hz, 10 Amp electrical service. Check the I.D. plate behind the front panel for voltage verification. The number of circuits varies depending on the model. All of the electrical connections are terminated at the field wiring terminal box for hook-up by a qualified electrician. We recommend each circuit be a dedicated branch circuit. However, if wired together the maximum load allowed is the sum of individual outlets plus the rating of the unit (i.e. 2 Amps remote, 8 Amps integral blower) not to exceed the 10A or 20A electrical service. The single point internal junction box is used for the connection of the lights, blower, and duplex outlets. Refer to the wiring diagram for your Protector PVC in *Appendix C: Protector PVC Specifications*.

The fume hood is required to be grounded to the MAINS protective earthing ground for safe operation. Using a ring terminal sized for a 10-24 machine screw, connect the MAINS ground conductor to the grounding lug marked with the protective earthing symbol, . Only MAINS ground conductors should be connected to the protective earthing ground lug, no other conductors should be connected to this grounding lug. Using wire nuts connect the MAINS supply conductors to the fume hood supply wires. Insure that the wires are connected as per the appropriate wire color codes for the input voltage. For 115V Phase (Hot) is black and Neutral is white, for 230V Phase1 is brown and Phase2 is blue. Refer to the wiring diagram for your Protector PVC in *Appendix C: Protector PVC Specifications*.

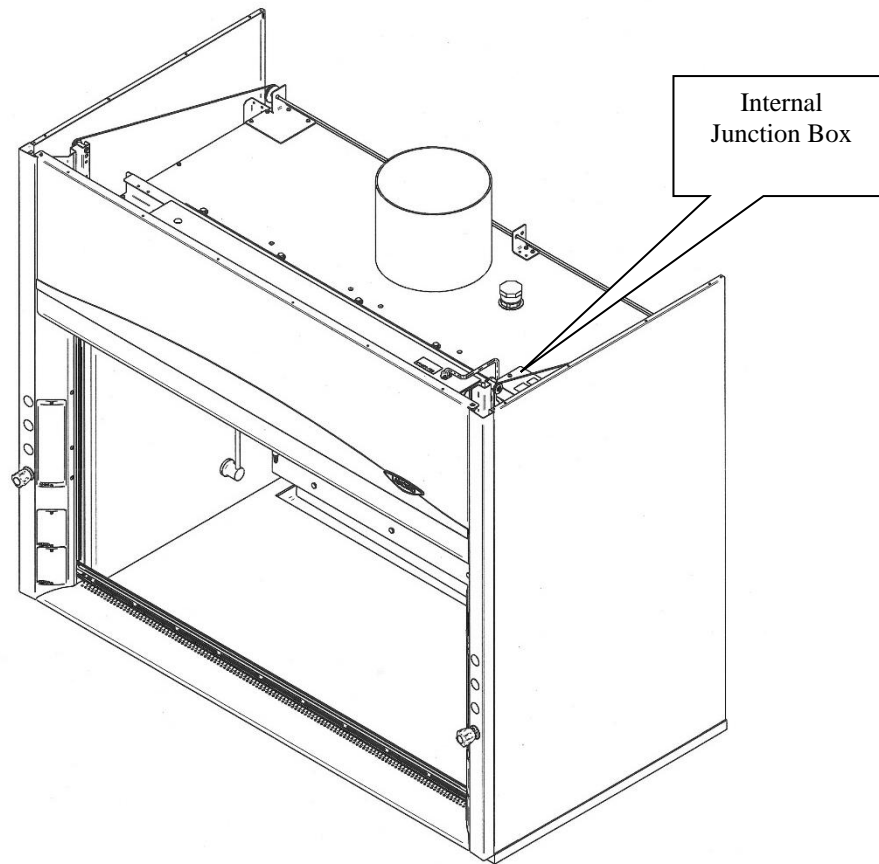
Figure 3-3

All wiring for the fume hood SHOULD be performed by a licensed electrician and conform to all local codes. In most cases, the hood will require the use of shielded conduit to protect the wiring into the hood. The grounding connection shall not be made to the terminal box cover.

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The LED light has been mounted outside the top liner panel and is sealed from vapors inside the hood structure. To change the fluorescent light bulbs in your hood, you must first remove the front panel from the hood. Next remove the knock out plugs holding the light fixture in place. Lift fixture up and replace any defective bulbs. Reverse order to reassemble.

Connecting the Service Lines to the Protector Fume Hood

The hoods with service fixtures have been plumbed from the valve to the hose connector or gooseneck for your installation convenience. Supply tubing shall be provided by the qualified installer. Tubing can enter the hood from above, through the back, or through the work surface to make these connections to the service fixtures.

NOTE: Inspect all fittings for leakage. Tighten the fittings slightly if needed.

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CAUTION: Do not use oxygen with any standard service fixture. Contact Labconco Customer Service for oxygen fixture information.

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Should access to the hood plumbing fixture bodies be required, remove the service access plate on the hood front corner posts by loosening their individual screws (see item 11, Figure A-1 in *Appendix A*). The valve body will now be fully exposed for any service work that may be necessary. The service fixtures supplied on your laboratory hood are designed for use with the following services:

- Air
- Hot Water
- Vacuum
- Cold Water
- Natural Gas – See caution below

WARNING: Contact Labconco Customer Service directly before using any service other than those listed above in these valves to assure full compatibility.

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CAUTION: Natural gas should be used only in the service fixture that has been pre-plumbed with brass tubing. Sulfur content of the gas could cause deterioration of standard copper supply lines.

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Sealing the Protector Hood to the Work Surface

When the hood has been set in place, ducted, wired, and plumbed, it should be sealed to the hardboard work top. Materials such as silicone sealants are recommended to seal the hood structure.

Certifying the Protector PVC Fume Hood

The combination of your laboratory hood, exhaust ductwork, and exhaust blower gives you the flexibility to change the airflow at the sash opening of your hood. To determine the actual face velocity at the sash opening, airflow velocity readings will need to be taken. This should be done across the sash opening of the hood in accordance with the *Industrial Ventilation Manual* section on laboratory hoods (see *Appendix E: References*). Labconco recommends an average face velocity at the sash opening of 100 feet per minute. Consult Chapter 2 for proper airflow volumes for your particular model.

Your Protector Fume Hood has been tested at the factory per ASHRAE 110-1995. All hoods achieve an “as manufactured rating” of less than 0.05 part per million (ppm) at 4 liters per minute (lpm); AM<0.05 (consult Labconco for individual fume hood ratings). For “field use” ASHRAE testing contact Labconco Sales Engineering Team or Customer Service for a certified on-site contractor.

NOTE: Face velocity profiles and smoke testing should be done periodically to ensure safe performance.

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AND SAFETY PRECAUTIONS

Performance Features

Your Protector PVC Perchloric Acid Fume Hood has been engineered to efficiently ventilate and remove Perchloric acid fumes. The Protector PVC Acid Digestion Fume Hood has been engineered to efficiently ventilate and remove acid fumes of most kinds (excluding Perchloric acid) including hydrofluoric acid (HF). The Protector PVC Acid Digestion Hood must not be used with Perchloric acid as it can react with the Lexan sash. The Protector PVC Fume Hoods are designed to provide a flow of air sufficient to extract the acid fumes from within the work area and keep them away from the operator to avoid ingestion, inhalation, or skin contact.

The hood interior is constructed of seamless PVC with coved interior corners. The hood also features an integral washdown system, which helps prevent the accumulation of perchlorates or acids. The smooth interior finish compliments the washdown system as there are no screws or hollow fasteners present that could harbor undesirable residue.

The Protector PVC Fume Hood has been engineered to allow you, the customer, the maximum work area and personnel safety. The by-pass air configuration on the Protector PVC Fume Hoods is totally dependent on air being supplied to the hood from its surrounding environment for proper operation.

As stated previously, the Protector PVC Perchloric Acid Hood is designed to handle Perchloric acid operations of all types, while the PVC Acid Digestion Hood is designed to handle acids of all kinds (excluding Perchloric acid) including hydrofluoric acid (HF). Usage of materials other than the specific acids mentioned above is not recommended. Perchloric acid is a strong oxidizing agent, and should not be used in combination with any organic material as the possibility of a hazardous chemical reaction or explosion could develop.

The integral washdown system in either PVC Fume Hood allows the operator the ability to wash behind the baffle and ensure that a build-up of Perchloric salts or other acids does not develop in that location. All of the washdown water from the internal spray nozzles is contained behind the baffle and does not interfere with normal operations being performed inside the hood. The washdown water is collected in the hood drain trough.

The hood liner **must** be washed down manually by the operator for **Perchloric acid**. The internal washdown system can be connected to the wash rings used in your hood exhaust system. The water used in the spray wash rings is collected in the hood drain trough once it has been allowed to flow down and cleanse the duct thoroughly.

WARNING: It is required that the PVC Perchloric Acid Hood be washed down thoroughly after every usage for 10 minutes, at least once a day.

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1. **Unique sash provides maximum visibility of 37.5" high while conserving energy by limiting sash travel to 28"**. Vertical-rising sash may be raised from a closed to 28" operating height. Exhaust volume and blower sizing are based on the 28" height. Optional sash stops are available to limit sash height and reduce energy usage. The PVC Perchloric Acid Hood has a glass sash specifically used for Perchloric Acid. The PVC Acid Digestion Hood has a Lexan sash specifically used for all acids (excluding Perchloric acid) including hydrofluoric acid (HF). **NOTE:** Hydrofluoric acid will etch or fog standard plate glass, but not Lexan plastic.
2. **By-pass airflow design** ensures containment as the sash moves.
3. **Large usable interior work depth and interior height of 48"** provides ample working space.
4. **Baffle** directs airflow to the rear of the interior to provide efficient airflow. The baffle may be removed for cleaning purposes only.
5. **Exterior access cover plates** are removable for easy access to plumbing valves when access through the sides is not available.
6. **Lift-Away™ front panel** provides easy access to electrical wiring, sash weights, and lighting fixtures.
7. **Energy efficient LED lighting** is located behind a laminated safety glass shield mounted to the top of the hood. The factory-wired lighting is serviceable from outside the hood cavity. Additionally, the long lasting 50,000 hour direct wired LED T8 bulbs are more energy efficient, utilizing approximately ½ the power of fluorescent bulbs. See Appendix A for wattage.
8. **Low-mounted, factory-wired light and blower switches** are ADA compliant.

Chapter 4: Performance Features & Safety Precautions

9. **Clean-Sweep™ Eco Air Foil** allows air to sweep the work surface for maximum containment. If needed, see Chapter 7 and Appendix A for the new PVC Eco Foil with PVC material that resists corrosion from mineral acids.
10. **Streamlined corner posts** provide maximum visibility and the flexibility to add services; services are most easily added at the factory, because of the PVC liner.
11. **All hoods are factory-prepared for up to 7 service fixtures.** Additional fixtures are available only as a factory special. Liner hole drilling is required to add another fixture in the field.
12. **One duplex electrical receptacle** is mounted on the right corner post. Receptacles are factory-wired to hood single point junction box. Additional electrical receptacles are available as factory or field installed.
13. **PVC interior liner** is seamless PVC with coved interior corners and integral work surface.
14. **Accessory Guardian™ Digital Airflow Monitor or Guardian Jr. Monitor** continuously monitors face velocity. An audio/visual alarm alerts the user to low airflow conditions. The right corner post is factory prepared to accommodate the Guardian Monitor (sold separately).
15. **Optional Energy-Reducing A-Style Combination Sash Models.** These combination sashes allow the operator to use the hood with sashes either half open vertically or horizontally to conserve energy. Optional sash stops prevent raising the vertical sash above the half-open and fully closed positions unless manually defeated by the operator.
16. **Outside frame of epoxy-coated steel and galvanized steel** is durable and corrosion resistant.
17. **Exhaust connection.** The hood features 12" (12.75" OD pipe) PVC exhaust connections sized to allow for a minimum static pressure loss through the hood structure while providing a good transport velocity through the exhaust system.
18. **Solid hardboard support** provided to support the integral PVC bottom.
19. **Optional Ceiling Enclosure Kits** are available for a decorative facade between the hood and the ceiling.
20. **Optional Sash Stops** provide a means of controlling the operating height of the sash.
21. **Washdown control fixture.** Located on the left hand side of the hood superstructure, this valve controls the water flow to the integral washdown spray nozzles. The nozzles are located behind the hood baffle and washdown areas in the hood, which are inaccessible without removing the baffle. The control fixture can also be set up to control the washdown system used in your exhaust ductwork. To do this, connect the water line from your external washdown system to the plugged connection provided on the top of the internal washdown assembly. For additional washdown valves and sizing, See Chapters 2 & 7.

22. **Acid Cabinet Vent Connection.** As a convenience, the PVC Hood has a 2" plugged connection on top of the hood to vent the acid cabinet below the fume hood. The acid vent kit is listed in Chapter 7 and ordered separately as part number 3591101.

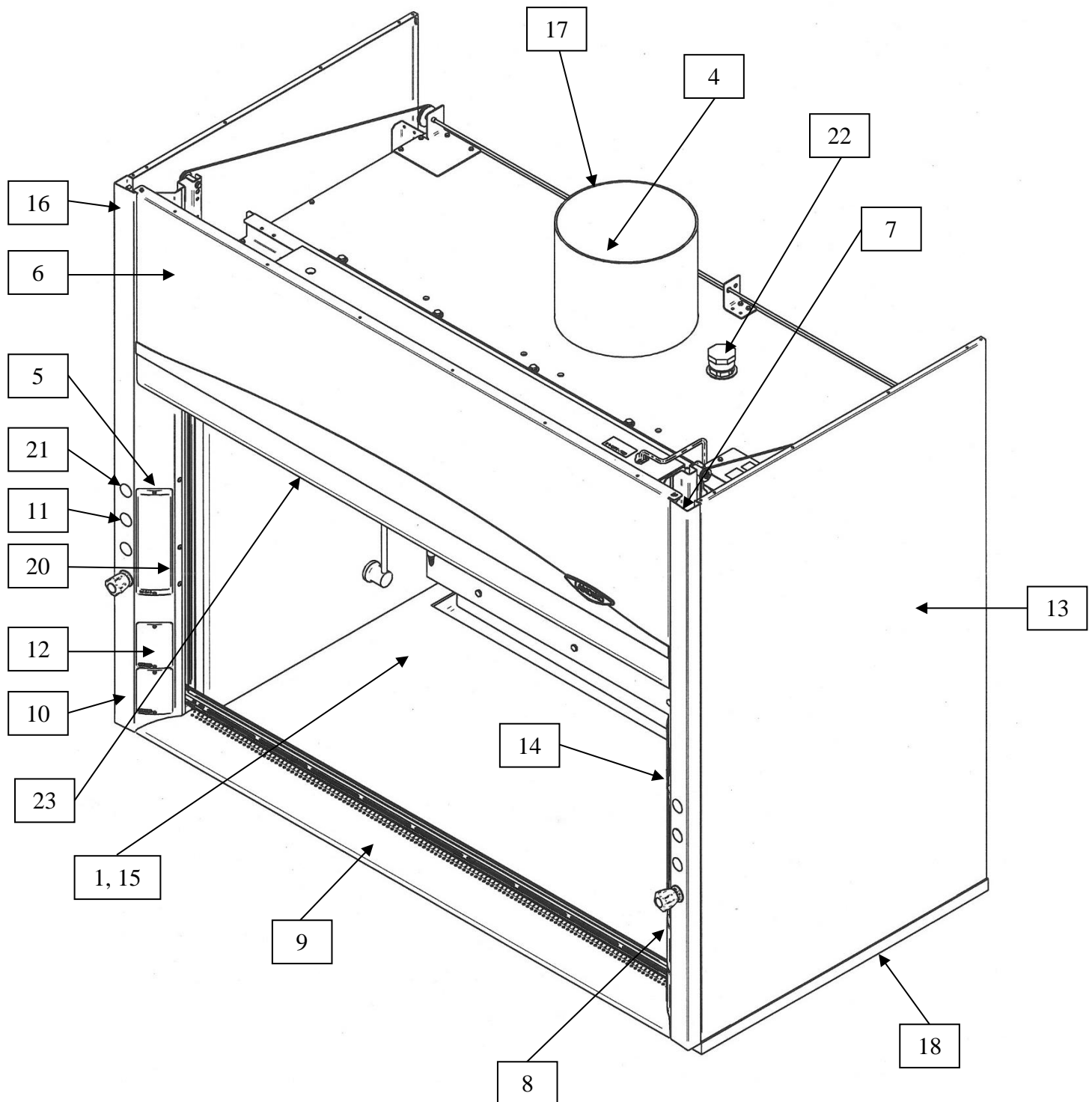
23. **Built-in Upper Bypass.** The PVC Hood includes an upper bypass block which proves to be useful when working with heat in digestions.

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26.  **CAUTION** – Flammable Gas.
AVERTIR – Gaz Inflammable.

Chapter 4: Performance Features & Safety Precautions**Figure 4-1**

Safety Precautions

Although the laboratory hood has been engineered to maintain optimum operator safety, caution should always be used while working in the hood. Prior to using the hood, check to make sure that the exhaust blower is operating and that air is entering the hood at its specified face velocity.

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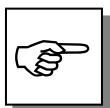
Do not place any hot items directly on the PVC Hood integral work surface.

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Always keep the ventilation system in operation when using any heat generating source inside the PVC Hood. Severe damage to the PVC Hood liner and PVC baffle can occur if the ventilation system is turned off before the heat load inside has been reduced to a safe level.

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USE GOOD HOUSEKEEPING IN THE HOOD AT ALL TIMES. CLEAN UP SPILLS IMMEDIATELY WITH A MILD DETERGENT. PERIODICALLY CLEAN HOOD INTERIOR, INCLUDING FLUORESCENT LIGHT GLASS PANEL. REPLACE BURNED OUT LIGHT BULBS TO MAINTAIN MAXIMUM ILLUMINATION.

DO NOT OVERLOAD THE WORK SURFACE WITH APPARATUS OR WORK MATERIAL. THE SAFE OPERATION OF THE LABORATORY HOOD IS BASED UPON HAVING PROPER AIRFLOW THROUGH THE STRUCTURE. DO NOT PLACE LARGE, BULKY OBJECTS SUCH AS BLOCK HEATERS, DIRECTLY ON THE HOOD

WORK SURFACE. INSTEAD, ELEVATE THE OBJECT 2" TO 3" ON BLOCKS TO ALLOW A FLOW OF AIR UNDER THE OBJECT AND INTO THE LOWER REAR BAFFLE EXHAUST SLOT. ENSURE BLOCKS ARE LEVEL AND SECURED IN PLACE.

Chapter 4: Performance Features & Safety Precautions

Blocking the bottom of the baffle at the rear of hood will change the airflow pattern in the hood causing turbulence and possible leakage at the face of the hood. (Don't store containers or supplies against baffles, as this will affect airflow through the hood).

Avoid placing your head inside hood. Keep hands out of hood as much as possible.

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Always work as far back in hood as possible. It is best to keep all chemicals and apparatus 6" inside the front of the hood.

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Perchloric acid use in PVC Perchloric Acid Hoods is recommended. Perchloric acid is not allowed in PVC Acid Digestion Hoods as it reacts with polycarbonate sashes.

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For Perchloric acid, always wash the hood down thoroughly after each usage or at least once a day for a period of 10 minutes.

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Safety requires regular washdown of fume hood interior, ductwork, and the movement of air preferably after each experiment. The baffle on this hood may be removed for access to all surfaces as a follow-up to washdown procedures. All deposits should be flushed away. Utilize lowest quantities of Perchloric acid or acids to fit procedural requirements.

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The use of Perchloric acid creates extreme inherent hazard because of its certain characteristics. Since there is a constant danger of explosion when using Perchloric acid, it is recommended that laboratories employing its use in any quantity become well informed of its characteristics. Only personnel fully cognizant with the properties of Perchloric acid and the hazards associated with it should perform Perchloric acid procedures.

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The PVC Perchloric Acid Hood is specially designed for use with Perchloric acid. It should not be used as a general-purpose laboratory fume hood. Do not work with organic materials in the PVC Perchloric Acid Fume Hood. Perchloric acid when used with organic material can cause a hazardous chemical reaction or explosion. Because of extreme hazards, miscellaneous work should not be performed in this hood. Do not store chemicals in a fume hood.

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Chapter 4: Performance Features & Safety Precautions

All apparatus used within the hood interior should have inorganic coatings and lubricants.

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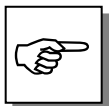
Although, the Protector PVC Perchloric Acid Fume Hoods conform to all generally accepted standards of design for this type of equipment, we accept no responsibility or liability for accidents that may possibly occur in the use of Perchloric acid in our hoods.

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Radioisotope materials are not recommended for use in PVC Hoods.

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AVOID CROSS DRAFTS AND LIMIT TRAFFIC IN FRONT OF THE HOOD. AIR DISTURBANCES CREATED MAY DRAW FUMES OUT OF THE HOOD.

The use of heat-generating equipment in this hood without the exhaust system operating properly can cause damage to the hood.

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The Protector PVC Laboratory Hoods should be certified by a qualified certification technician before initially used. The hood should be re-certified whenever it is relocated, serviced or at least annually thereafter.

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Ensure that the hood is connected to electrical service in accordance with local and national electrical codes. Failure to do so may create a fire or electrical hazard. Do not remove or service any electrical components without first disconnecting the hood from electrical service. Proper operation of the fume hood depends largely upon the hood's location and the operator's work habits. Consult the *Reference Manual* in *Appendix D*.

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If the unit is not operated as specified in this manual it may impair the protection provided by the unit.

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Do not position the fume hood so that it is difficult to operate the main disconnect device.

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USING THE PROTECTOR PVC

S

Operating the Vertical-Rising Glass Sash

Because of the Protector PVC Perchloric Acid Hood's counterbalanced sash mechanism, it will take only a few pounds of force to move the sash up or down, and you can operate the sash smoothly with one or two hands positioned anywhere along the handle. The glass sash is suitable for use with Perchloric acid. The vertical-rising sash may be raised to a maximum 28" operating height. The airflow requirements should be sized for the 28" operating height; if using sash stops then the airflow requirements can be reduced by approximately 40% at 18" or approximately 50% at 15".

S

Operating the Vertical-Rising Lexan Sash

Because of the Protector PVC Acid Digestion Hood counterbalanced sash mechanism, it will take only a few pounds of force to move the sash up or down, and you can operate the sash smoothly with one or two hands positioned anywhere along the handle. The Lexan sash prevents fogging caused by hydrofluoric acid (HF). The Lexan sash is suitable for acids other than Perchloric acid, which can react with the Lexan plastic. The airflow requirements should be sized for the 28" operating height, if using sash stops then the airflow requirements can be reduced by approximately 40% at 18" or approximately 50% at 15".

S

Operating the Blower

Your Protector PVC Fume Hood utilizes a remote style blower, which can be activated by turning the blower switch to "ON." You can validate the hood performance by watching smoke drawn into the hood face opening.

S

Operating the Lights

Your Protector PVC Fume Hood utilizes a factory-wired LED light to illuminate the hood interior (see Appendix A). Simply turn the light switch to "ON" to operate.

S

Operating the Washdown Control Valve

Your Protector PVC Fume Hood has a washdown control valve located on the upper left-hand side that controls water to the washdown spray nozzles. The nozzles are located behind the hood baffle and will washdown areas in the hood, which are inaccessible without removing the baffle. The washdown control system can also be set up to control the washdown rings used in your exhaust ductwork. This is done by connecting the water line to the plugged connection on top of the washdown hood system or ordering additional washdown valves found in Chapter 2 and Chapter 7.

Working in Your Protector PVC Fume Hood

Planning

- Thoroughly understand procedures and equipment required before beginning work.
- Arrange for minimal disruptions, such as room traffic or entry into the room while the hood is in use.

Start-up

Labconco follows the recommendations of the American National Standard Institute with regards to fume hood exhaust systems. A full list of requirements can be found in the ANSI Standard No. Z9.5 *Laboratory Ventilation*. If you are unsure of the proper use of the exhaust blower for your fume hood, contact your safety officer or industrial hygienist before making any modifications to the blower's operational status (on/off/low, etc.). Keeping the exhaust system under constant negative pressure is an excellent way to ensure hazardous fumes in the fume hood do not enter the laboratory space inadvertently.

- Turn on the light.
- Slowly raise the sash and verify the fume hood has proper airflow.
- Check the baffle air slots for obstructions.
- Allow the hood to operate unobstructed for two minutes.
- Wear a long sleeved lab coat and rubber gloves. Use protective eyewear. Wear a protective mask if appropriate.

Loading Materials and Equipment

- Only load the PVC Hood materials required for the procedure. Do not overload the hood. Only use Perchloric acid in the PVC Perchloric Acid Fume Hood.
- Radioisotope materials are not recommended for use in these PVC Hoods. Consult your Safety Officer should you have any questions.
- Do not obstruct the front air foil (sill), or rear baffle slots.
- Large objects should not be placed close together and spaced above the liner bottom to permit airflow to sweep under the equipment.
- After loading the hood, wait one minute to purge airborne contaminants from the work area.

Chapter 5: Using the Protector PVC

Work Techniques

- Keep all materials at least 6 inches inside of the sash and at least two inches from the side walls and perform all contaminated operations as far to the rear of the work area as possible.
- Keep all clean and contaminated materials in the work area separate for clear identification.
- Avoid using techniques or procedures that disrupt the airflow patterns of the hood.

Final Purging

- Upon completion of the work, the hood should be allowed to operate for two minutes undisturbed, to purge airborne contaminants from the work area before closing the sash.

Unloading Materials and Equipment

- Objects in contact with contaminated material should be surface decontaminated before removal from the hood.
- All open trays or containers should be covered before being removed from the hood.

Final Washdown Procedure

- Water washdown or flooding the surfaces with water is standard practice and results in reduction of contamination on the PVC liner material due to Perchloric acid salts or acids.
- Proper operation of your PVC Fume Hood requires that you wash the fume removal system down sufficiently to ensure that there is no build up of Perchloric acid salts or acids within the entire system.
- Normal washdown procedures suggest that you run the wash system in the exhaust duct with the fan on for approximately 1-2 minutes, and then with the fan off for approximately 10 minutes to thoroughly clean this portion of the exhaust system.
- In addition to this, the internal hood spray nozzles should be activated for approximately 10 minutes after the completion of your procedures to clean behind the baffle in the hood. Work areas in front of the baffle in the hood will need to be manually cleaned by the operator. Because the volume of Perchloric acid salts or acids can vary greatly due to specific procedures and workloads, you should consult your in-house Safety Officer for specific rules and frequency requirements for washing down your total system.

Shutdown

- Only turn off the light and close the sash if final washdown is complete.

CHAPTER 6

MAINTAINING THE PROTECTOR PVC

Now that you have an understanding of how to work in the fume hood, we will review the suggested maintenance schedule and the common service operations necessary to maintain your fume hood for peak performance.

Service Safety Precautions

- If performing any electrical maintenance, always disconnect the power at the main disconnect.
- If performing decontamination inside the fume hood, consult your safety officer for proper personal protective equipment and procedure.
- Since some service operations require a step ladder, always use proper safety and consult your safety officer
- If performing maintenance on any service lines, always shut off the supply first.
- Some removable components may be heavy, follow safe-lifting guidelines.
- Verify all components are installed correctly with performance verified before conducting normal operations.

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Chapter 6: Maintaining the Protector PVC

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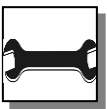
Weekly

- Using ordinary dish soap to clean the surface inside of the fume hood, and the work surface.
- Using an appropriate glass cleaner, clean the sash and all glass surfaces.
- Operate the fume hood blower, noting the airflow velocity through the hood using a source of visible smoke.

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Biannually

- The sash assembly should be checked to ensure proper operation and to make sure there are no signs of abnormal wear on the sash pulleys, cables and clamps.

Routine Service Operations

Front Panel Removal:

1. Simply lift the front panel up and then away from the hood to provide access to the top.

Changing the Lamp:

1. Turn light switch to “OFF”.
2. Remove the front panel as noted earlier.
3. Reach over the front header of the hood and remove knock out plugs at both ends of fixture. Lift fixture up.
4. Remove the lamp by pushing it out of the spring-loaded lamp socket and swinging it out of the other lamp socket.
5. Install the new lamp by reversing the removal procedure.

CHAPTER 7

MODIFYING THE PROTECTOR PVC

There are several ways to modify the fume hood for your individual requirements. These include the addition of work tops, service fixtures, air monitor, electrical duplex outlets, ceiling enclosures, and rear panels.

Installing Hardboard Support Surface

Your Protector PVC Fume Hood requires a hardboard support surface to properly support the integral work surface of the one-piece PVC liner. The hardboard support is shipped and included with the PVC Fume Hood. Be sure to secure the hardboard support per the instructions in Chapter 3.

Installing PVC Eco Foil Made from PVC Material

Your Protector PVC Fume Hood normally uses a painted stainless steel lower Eco Air Foil. However, some customers require a PVC Eco Air Foil manufactured from PVC material. If a PVC Eco Air Foil is needed to enhance corrosion resistance from acids, then please see Appendix A, items 13E through 13N to order. Order your size from 13E to 13H and all hardware 13I through 13N.

Installing Ceiling Enclosures Above the Fume Hood

Your Protector Fume Hood has mounting holes to accept a ceiling enclosure to close off the area between the top of the hood and the ceiling. Contact Labconco Customer Service for ordering information. Labconco offers both non-adjustable fixed height ceiling enclosures as well as adjustable height ceiling enclosures per the charts below. Other special sizes are available upon request.

3-Sided Ceiling Enclosure Kits

The panels extend above the top of the hood to the ceiling to hide exposed ductwork, plumbing and wiring. For any 3 ft to 8 ft Protector XL Benchtop, Premier, XStream, Stainless Steel or PVC Fume Hood. Each kit includes three glacier white powder-coated steel ceiling enclosure panels, one for the front of the hood and one for each side of the hood.

Configure Your Non-Adjustable Fixed Height Ceiling Enclosure Kit Catalog Number

Use this key to configure the seven digit catalog number to order your 3-Sided **Fixed Height** Ceiling Enclosure Kit. For example, **9550511** is a 3-Sided Ceiling Enclosure Kit for a 5' Protector PVC Hood with a fixed height of 12".

9

5

5

0

STEP 1. Select the **nominal width** of your fume hood. This number is the fifth digit of your catalog number.

3 = 3' (XL Benchtop only)

4 = 4'

5 = 5'

6 = 6'

7 = 7' (XL Benchtop only)

8 = 8'

STEP 2. Select the **depth** based on the model chosen. Choose 0 for Premier and 1 for XStream. Choose 1 for PVC. Choose 0, 1 or 2 for the correct depth of the XL Benchtop model. Choose 0 or 1 for the correct depth of Stainless Steel. This is the sixth digit of your catalog number.

No.	Ext. Depth	Model
0	31.7"	Premier, XL Benchtop, Stainless Steel
1	37.7"	XStream, XL Benchtop, XL Floor-Mounted, Stainless Steel, PVC
2	43.7"	XL Benchtop, XL Floor-Mounted

STEP 3. Choose the correct **height** that matches the number. This height should match the height between the top of the hood and the building ceiling. This is the seventh digit of your catalog number.

No.	Height
0	6"
1	12"
2	18"
3	24"
4	30"
5	36"

Chapter 7: Modifying the Protector PVC

Configure Your Adjustable Height Ceiling Enclosure Kit Catalog Number

Use this key to configure the seven digit catalog number to order your 3-Sided **Adjustable Height** Ceiling Enclosure Kit. For example, **9425511** is a 3-Sided Ceiling Enclosure Kit for a 5' Protector PVC Hood with a height range of 14-19".

9 **4** **2** **5**

STEP 1. Select the **nominal width** of your fume hood. This number is the fifth digit of your catalog number.

3 = 3' (XL Benchtop only)
4 = 4'
5 = 5'
6 = 6'
7 = 7' (XL Benchtop only)
8 = 8'

STEP 2. Select the **depth** based on the model chosen. Choose 0 for Premier and 1 for XStream. Choose 1 for PVC. Choose 0, 1 or 2 for the correct depth of the XL Benchtop model. Choose 0 or 1 for the correct depth of Stainless Steel. This is the sixth digit of your catalog number.

No.	Ext. Depth	Model
0	31.7"	Premier, XL Benchtop, Stainless Steel
1	37.7"	XStream, XL Benchtop, XL Floor-Mounted, Stainless Steel, PVC
2	43.7"	XL Benchtop, XL Floor-Mounted

STEP 3. Choose the correct **height range** that matches the number. This height range should match the height between the top of the hood and the building ceiling. This is the seventh digit of your catalog number.

No.	Height Range
0	11-14"
1	14-18.6"
2	18.6-24.4"

Note: The previous generation listed as 9414-W-D-H is presented below to aid product service.

9 **4** **1** **4** **W** **D** **H**


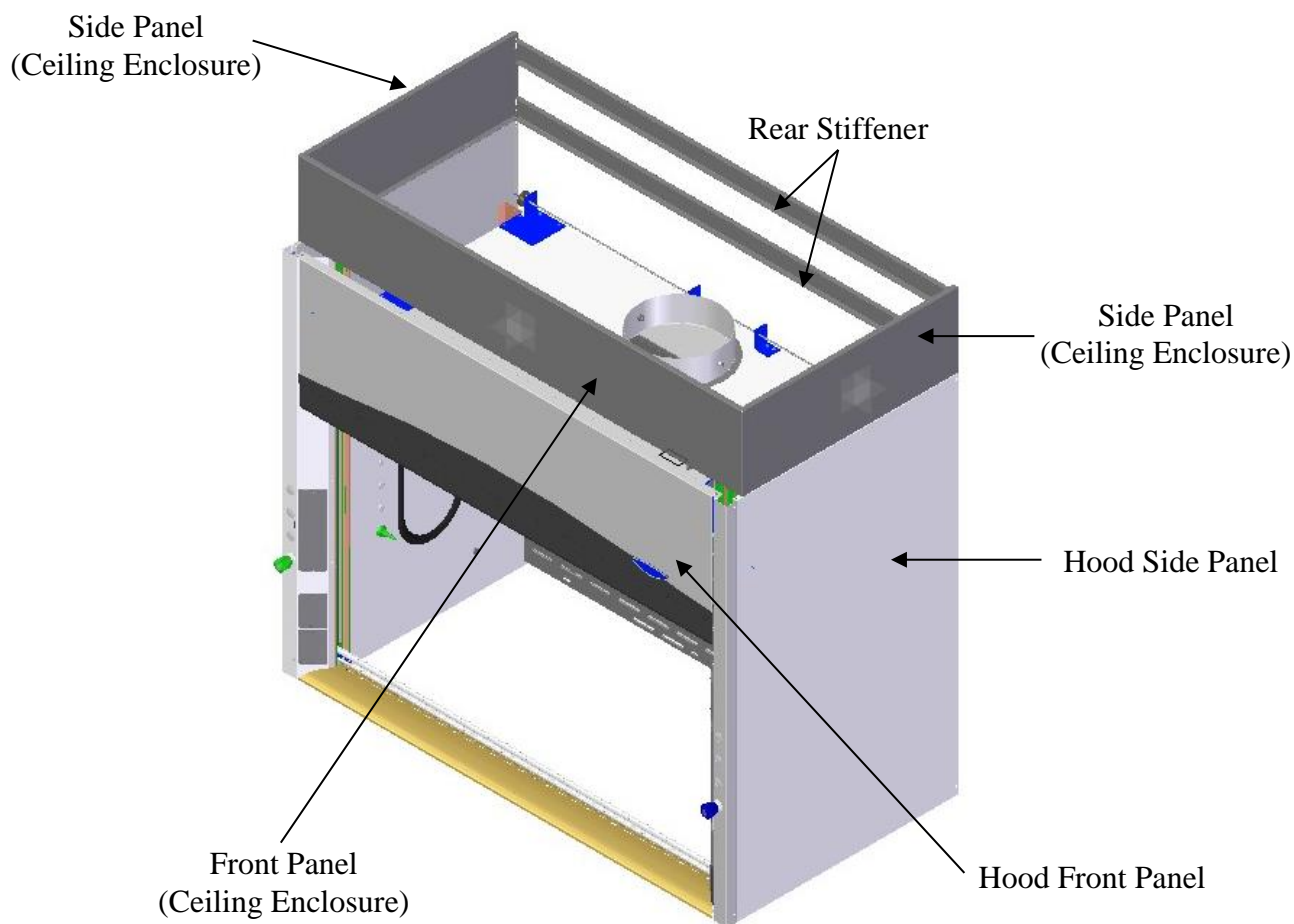
Figure 7-1

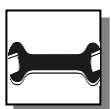
Installing Rear Panels Behind the Fume Hood

Your Protector Fume Hood can be modified to add a rear panel behind the fume hood when the fume hood is placed on an island (not available on Pass-Through Fume Hood). Contact Labconco Customer Service for ordering information.

Installing Guardian™ Digital Airflow Monitor or Guardian™ Airflow Monitor

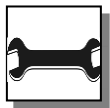The Guardian Airflow Monitors continuously monitor face velocity through the fume hood opening. The fume hood right corner post is factory prepared to mount either monitor. Contact Labconco Customer Service to order.

Sash Stop Kit – Field Installation (P/N 9410300)

The sash stop kit restricts how far a vertical-rising sash may be opened. This small plastic device may be easily field installed on the fixture corner post of any fume hood.

Installing an Electrical Duplex Outlet

Your Protector Fume Hood can be ordered with duplex outlets, however, if you ordered a model without an electrical duplex outlet you can have one installed in the field by a qualified electrician. Contact Labconco Customer Service for ordering information. (Not acceptable on explosion-proof hoods).

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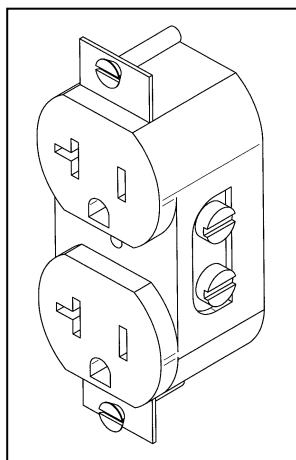
Figure 7-2

Installing the Acid Vent Kit from the Acid Cabinet

Order 3591101 or 3591102 Acid Vent Kit for PVC hoods to install from the Acid Cabinet below to the top of the PVC Hood. Remove the plugged 2" connection and route the tubing.

Installing Additional Washdown Valve

Order 9808310 Washdown Valve and follow the sizing guideline from Chapter 2.

CHAPTER 8

TROUBLESHOOTING

Refer to the following table if your fume hood fails to operate properly. If the suggested corrective actions do not solve your problem, contact Labconco for additional assistance.

PROBLEM	CAUSE	CORRECTIVE ACTION
Remote blower and lights won't operate	Wires not connected at junction boxes or switches.	Check connection of switches.
		Check connection to control box on top of unit.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Remote blower won't operate, but lights work	Blower wiring is disconnected.	Inspect blower wiring and switch.
	Belt broken.	Replace belt.
	Blower motor is defective.	Replace blower motor.
Fume hood blower operates but lights will not operate	Lamp not installed correctly.	Inspect lamp installation.
	Lamp is defective.	Replace lamp.
	Lamp circuit breaker in building is tripped.	Reset the lamp circuit breaker.

Chapter 8: Troubleshooting

PROBLEM	CAUSE	CORRECTIVE ACTION
Fume hood blower operates but lights will not operate	Lamp wiring is disconnected.	Inspect lamp wiring.
	Defective lamp ballasts.	Replace lamp ballasts.
Contaminants escape outside of fume hood	Improper user techniques for the fume hood.	See “Certifying the Hood” Chapter 3 and “Safety Precautions” Chapter 4 sections in the manual. (Ref. Appendix D).
	Restriction of the baffle air slots or – blockage of the exhaust outlet.	Remove baffles to ensure that all air slots and the exhaust outlet are unobstructed.
	External factors are disrupting the fume hood airflow patterns or acting as a source of contamination.	See “Location Requirements” Chapter 2, “Certifying the Hood” Chapter 3, and “Safety Precautions” Chapter 4 sections of this manual. (Ref. Appendix D)
	Fume hood has improper face velocity.	Have fume hood re-certified and check remote blower exhaust system. Hood should have average face velocity of 100 fpm.
Vertical sash no longer operates smoothly	Cable is frayed or plastic protection is damaged.	Inspect cable and replace cable if worn or damaged immediately; otherwise injury could result.
	Pulley bearing is damaged.	Replace pulley, bearing or add grease.
	Cable has slipped off the pulleys.	Re-install, cable must be replaced immediately if damaged.
	Weight has broken pulleys.	Replace weight pulleys.
Combination A-Style sash no longer operates smoothly	Horizontal-sliding glass panels have come off the tracks.	Re-install horizontal-sliding glass on tracks.

PROBLEM	CAUSE	CORRECTIVE ACTION
Electrical duplex outlets no longer have power	Vertical-rising sash frame is distorted.	Place horizontal sliding glass symmetrically and pull sash down to air foil. Straighten damaged frame.
	Cable is frayed or has slipped off the pulleys.	Re-install, cable must be replaced immediately if damaged.
	Wires not connected or faulty duplex.	Check wire connection or replace duplex.
	Circuit breakers tripped in building electrical supply.	Reset circuit breakers.
Service valves no longer operate	Faulty building supply.	Inspect building supply shut off valves and appropriate pressures below the recommended 40 PSI.
	Valve no longer operates.	Replace valve and check for leaks.
	Supply line or outlet line has leaks.	Inspect line for leaks and fix any leaking plumbing connections.
Spray Nozzles no longer work	Poor water supply	Inspect building water supply.
	Clogged nozzle	Replace spray nozzles.

APPENDIX A

PROTECTOR PVC COMPONENTS


Illustration A-1 indicate the location of the following service parts:

Protector PVC Replacement Parts

Item	Quantity	Part No.	Description
1A	1	9823700	Valve, Labconco (Water) 1/4" Compression Fitting
1B	1	9823701	Valve, Labconco (Water) 3/8" Compression Fitting
1C	1	9817000	Valve, Labconco 1/4" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1D	1	9817001	Valve, Labconco 3/8" Compression Fitting (AIR, GAS, VAC, NIT, etc.)
1E	1	9823702	Valve, Labconco Deionized 1/4" Compression Fitting
1F	1	9823703	Valve, Labconco Deionized 3/8" Compression Fitting
1G	1	9818000	Nut, Valve Mounting (Labconco)
2A	1	9826800	WaterSaver Valve/Gooseneck – GRN
2B	1	9826801	WaterSaver Valve/Connector (VAC) – YEL
2C	1	9826802	WaterSaver Valve/Connector (AIR) – ORG
2D	1	9826803	WaterSaver Valve/Connector (GAS) – BLU
2E	1	9826805	WaterSaver Valve/Connector (HOT WATER) – RED
2F	1	9826806	WaterSaver Valve/Connector (CW) – GRN
2G	1	9826807	WaterSaver Valve/Connector (STEAM) – BLK
2H	1	9826808	WaterSaver Valve/Connector (NITROGEN) – BRN
2I	1	9826809	WaterSaver Valve/Connector (OXYGEN) – LIGHT GREEN
2J	1	9826810	Swivel Gooseneck only – GRN
2K	1	9826812	Swivel Gooseneck only – WHITE
3	1	9818700 thru 08	Knobs (GRAY, GRN, BLU, ORG, YEL, RED, WHT, BLK, BRN)
4A	1	9818800	Hose Barb, GRAY (NEUTRAL OR ARGON) – NOT SHOWN
4B	1	9818801	Hose Barb, GREEN (COLD WATER) – NOT SHOWN
4C	1	9818802	Hose Barb, BLUE (GAS) – NOT SHOWN
4D	1	9818803	Hose Barb, ORANGE (AIR) – NOT SHOWN
4E	1	9818804	Hose Barb, YELLOW (VACUUM) – NOT SHOWN
4F	1	9818805	Hose Barb, RED (HOT WATER) – NOT SHOWN
4G	1	9818806	Hose Barb, WHITE (DEIONIZED WATER) – NOT SHOWN
4H	1	9818807	Hose Barb, BLACK (NEUTRAL OR STEAM) – NOT SHOWN
4I	1	9818808	Hose Barb, BROWN (NITROGEN) – NOT SHOWN
4J	1	9819000	Nut, Hose Barb – NOT SHOWN
5	1	9825500	Label, Knob (contains all the labels)
6	1	9818900	Lens, Knob
7A	1	9947100, 01, 02	115V Duplex Receptacle (GRAY) Right, Left 4' - 6', Left 8' w/ wires
7B	1	9818200	Cover Plate 115V Duplex
7C	1	9947103, 04, 05	115V GFCI Duplex Receptacle (GRAY) Right 4' - 6', Left 8' w/ wires
7D	1	9818100	Cover Plate, 115V GFCI
7E	1	9818300	Cover Plate, Blank
8A	1	1297503	LED, Direct Wired T8 x 3' – newest models use on 4' & 8' Hoods – 12 watts each
8B	1	1297504	LED, Direct Wired T8 x 4' – newest models use on 5' & 6' Hoods – 22 watts each
8C	1	9945300	LED Light Fixture, 36" – use on 4' Hoods
8D	1	9945301	LED Light Fixture, 48" – use on 5', 6' Hoods

Appendix A: Protector PVC Components

Item	Quantity	Part No.	Description
8E	1	9945302	LED Light Fixture, 36" Combo – use on 8' Hoods
10A	1	1302301	Switch, Rocker
10B	1	1327500	Switch, Plug (Fills cutout when switch is not used)
11	1	9818400	Access Cover
12A	1	9409800	Side Panel, 24" internal deep hoods
12B	4	1885308	Screw, Machine #10-24 x .50 Phillips
13A	1	9466400	Eco-Foil 4' (powder coated stainless p/n)
13B	1	9466401	Eco-Foil 5' (powder coated stainless p/n)
13C	1	9466402	Eco-Foil 6' (powder coated stainless p/n)
13D	1	9466403	Eco-Foil 8' (powder coated stainless p/n)
13E	1	9466420	Eco PVC Foil, 4' for PVC Hoods (optional new style made from PVC)
13F	1	9466421	Eco PVC Foil, 5' for PVC Hoods (optional new style made from PVC)
13G	1	9466422	Eco PVC Foil, 6' for PVC Hoods (optional new style made from PVC)
13H	1	9466423	Eco PVC Foil, 8' for PVC Hoods (optional new style made from PVC)
13I	4	9491901	Bracket, Airfoil Support – PVC Foil on PVC Hoods
13J	8	1937317	KEPS Nut #10-24
13K	8	1889308	Screw, Machine Phil Ph #10-24 x .50 SS
13L	4	1894712	Screw, Self-Drilling Hex #10 x .75
13M	4	1911210	Washer, Fender #10 SS
13N	4	1598500	Cover-Hinged Plastic
14	2	1850000	Pulley, Front, 2" Dia. Nylon
15A	2	4949902	Cable, Sash 130" – NOT SHOWN
15B	2	9414011	Cable Replacement Kit 30"
15C	1	9545800	Weight Support Bracket Kit
16	2	9709300	Plastic Pulley, (Rear 2")
17	2	9713300	Bumper, Rubber (lower sash bumper) – NOT SHOWN
18	4	1934100	Bronze Bearing, Flanged Rear – NOT SHOWN
19	4	1920100	Clamp, Cable Replacement – NOT SHOWN
20	2	9935800	Threaded Connector (to attach weight to cable) – NOT SHOWN
21	1	9946300 or 01	Wiring Harness, Main
22	1	9409601, 02, 03, 05	Front Panel, 4', 5', 6', 8'
23	1	1487100	Spray Nozzle PVC – NOT SHOWN
24	1	9817001	Washdown Control Valve
25	1	1549100	Plastic Tubing 3/8 OD x 1/4 ID (125 psi W.P.) (Rigid Polyethylene – PE) – NOT SHOWN

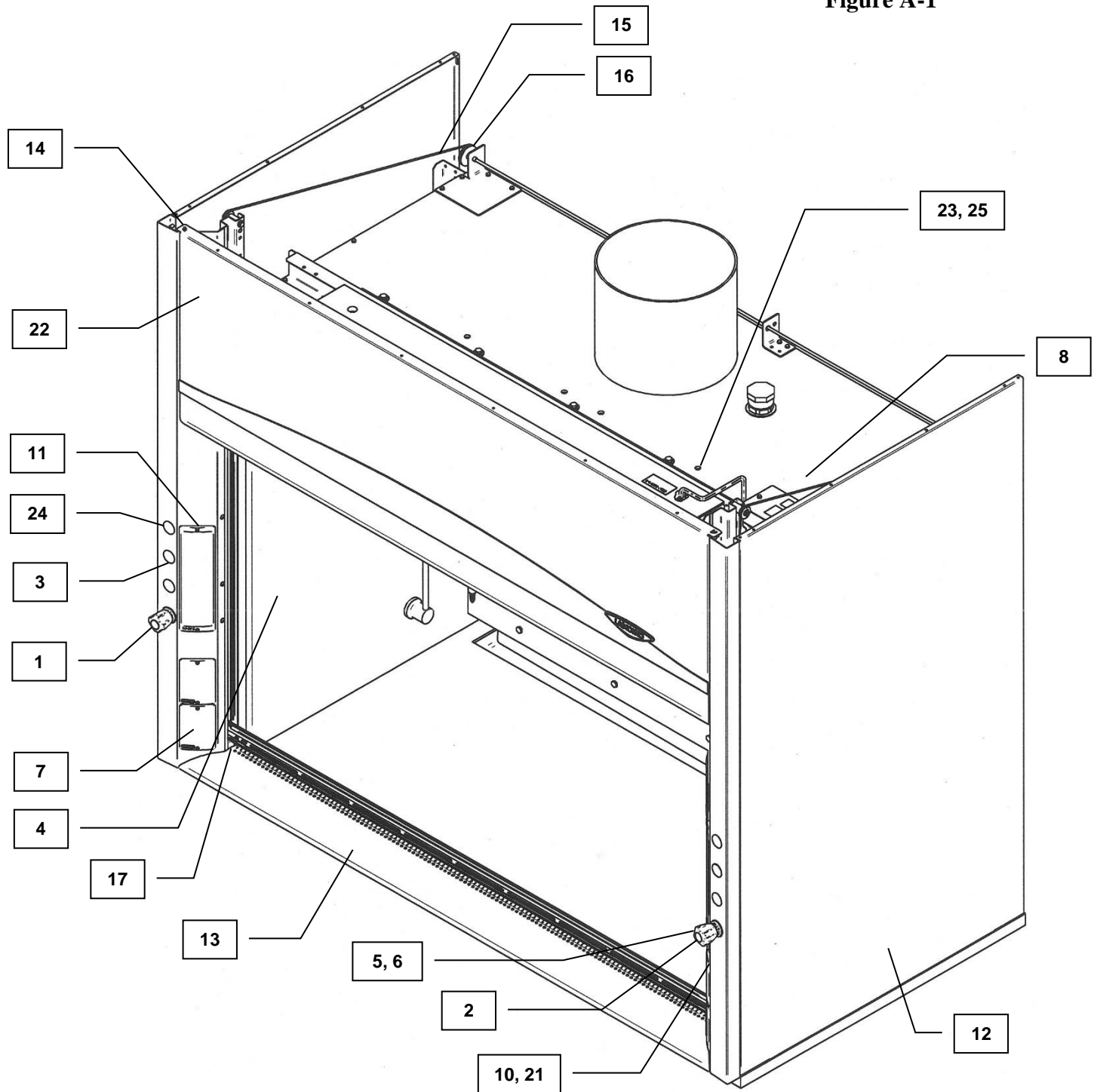Appendix A: Protector PVC Components**Figure A-1**

APPENDIX B

PROTECTOR PVC DIMENSIONS

PVC Model Widths

	4'	5'	6'	8'
A	48.00	60.00	72.00	96.00
B	38.25	50.25	62.25	86.25

Dimensions in inches. CAD layouts available upon request.

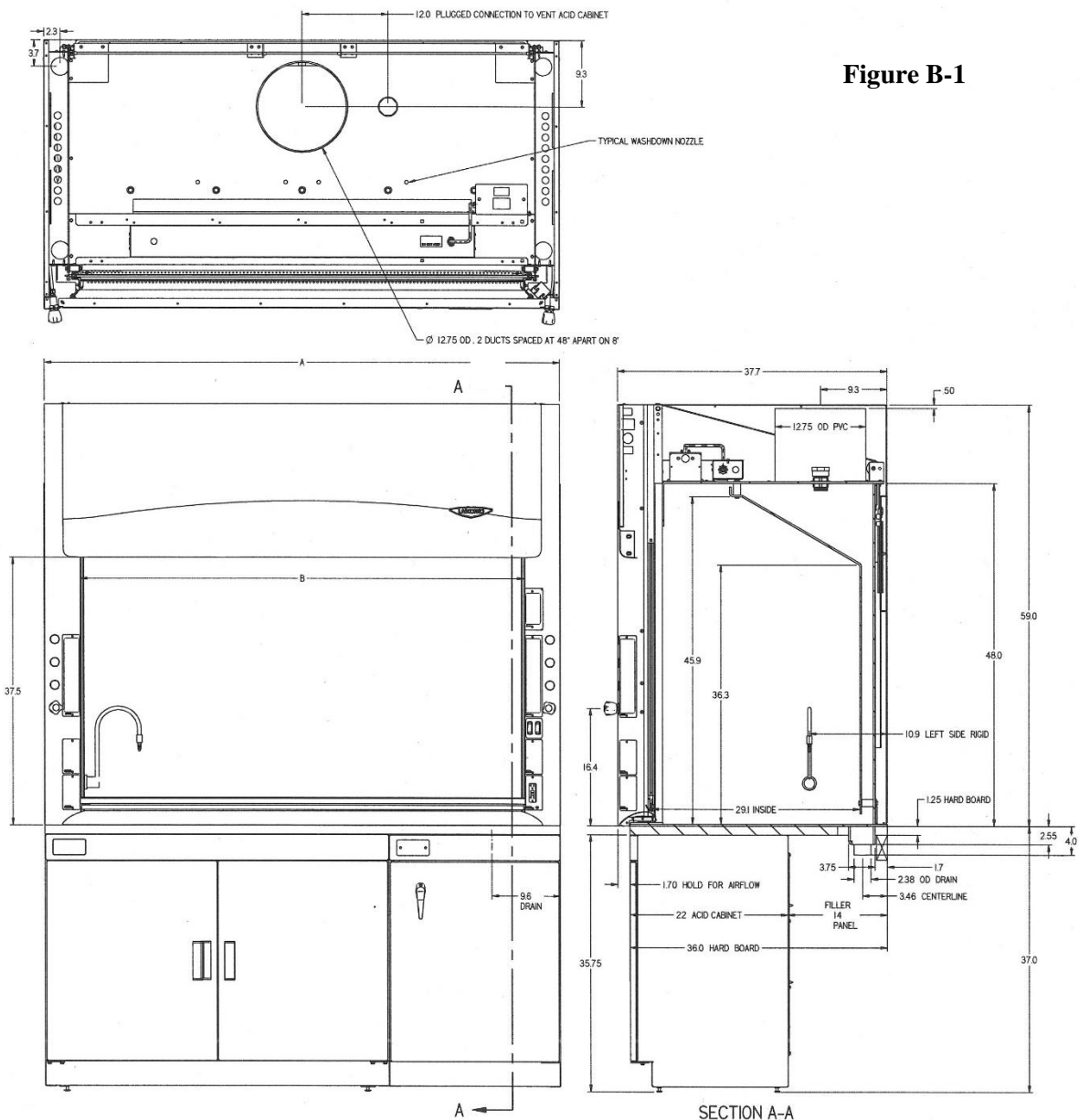
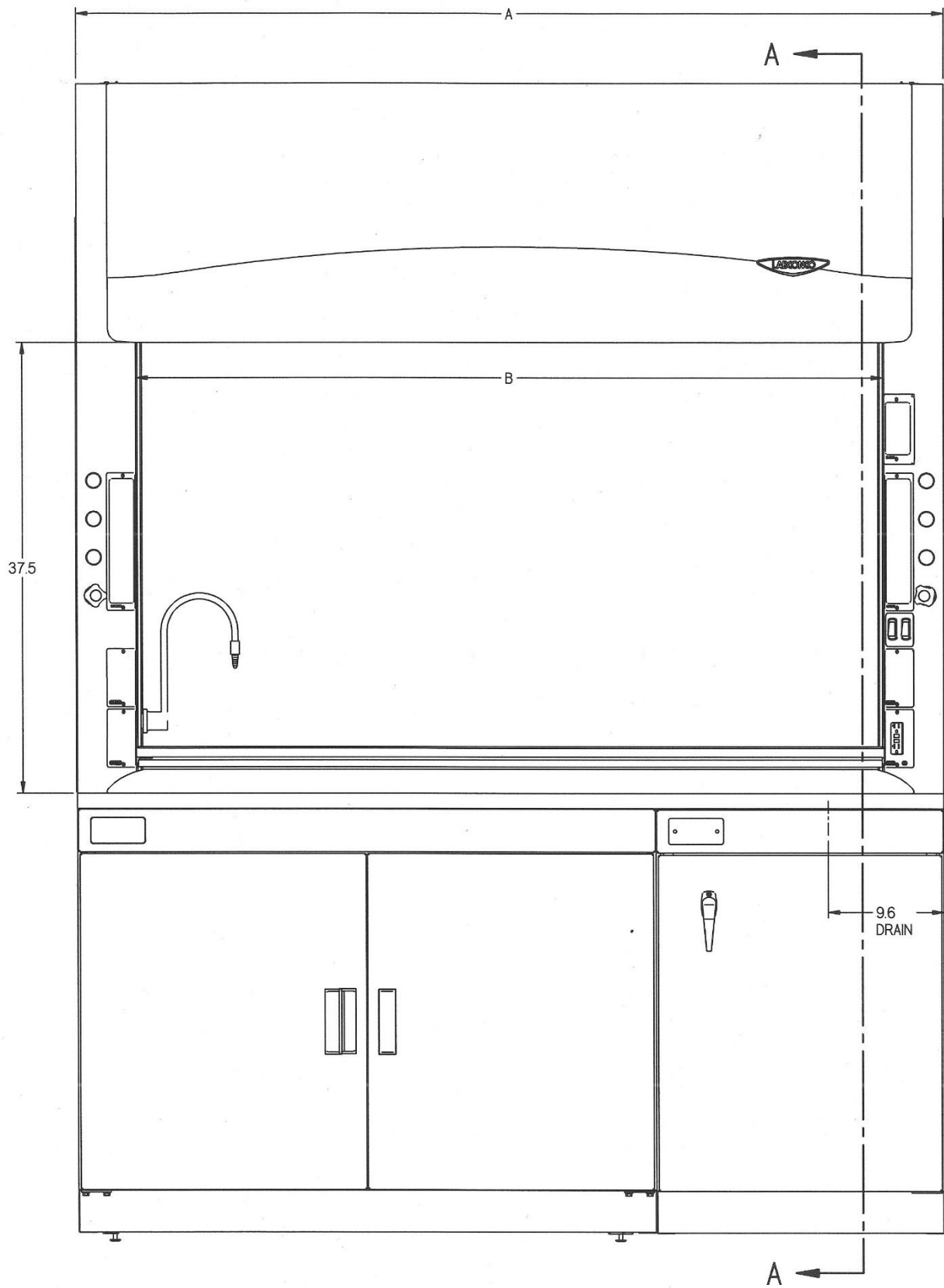
Figure B-1

Appendix B: Protector PVC Dimensions**Figure B-2**

Appendix B: Protector PVC Dimensions

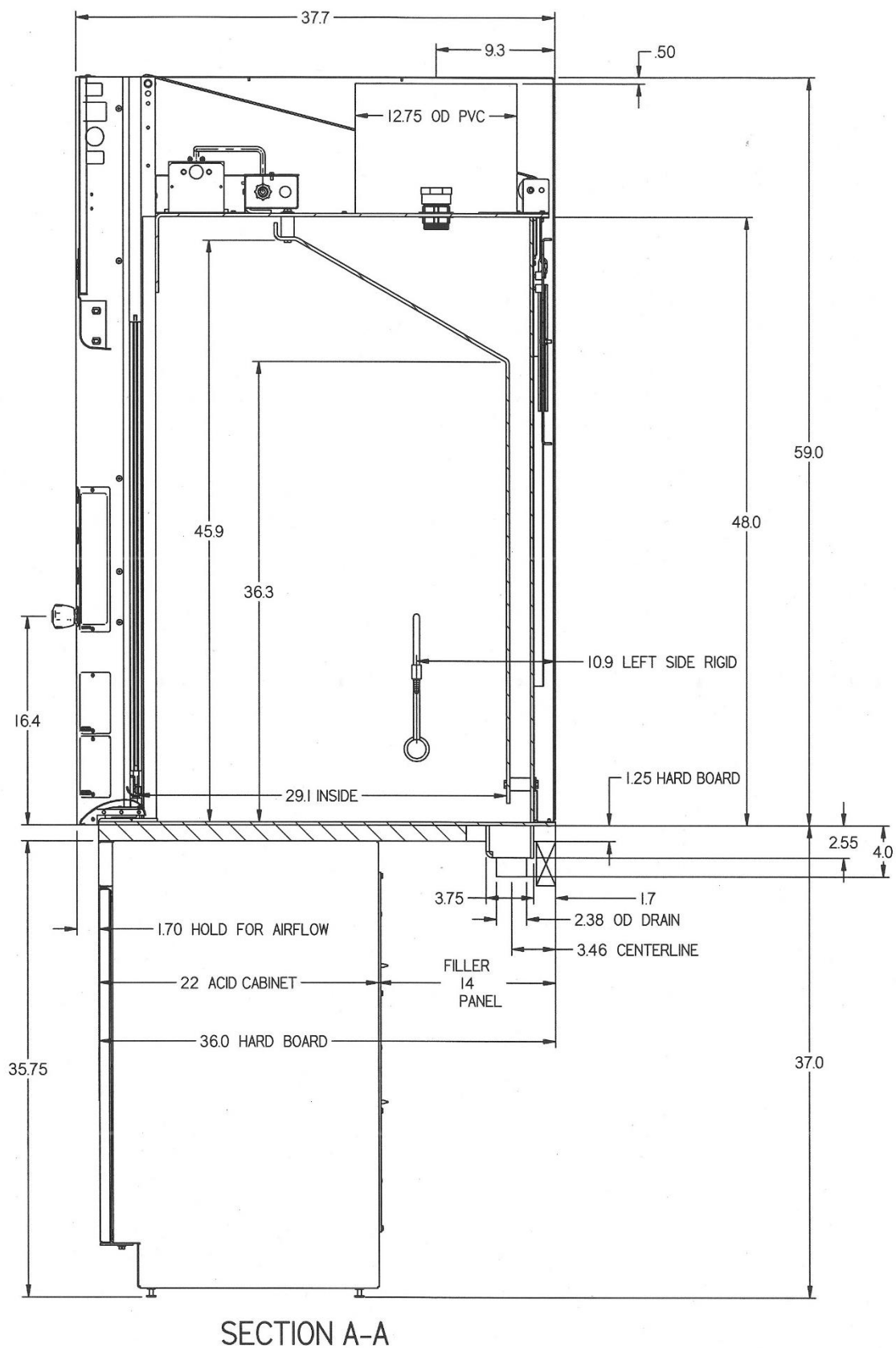
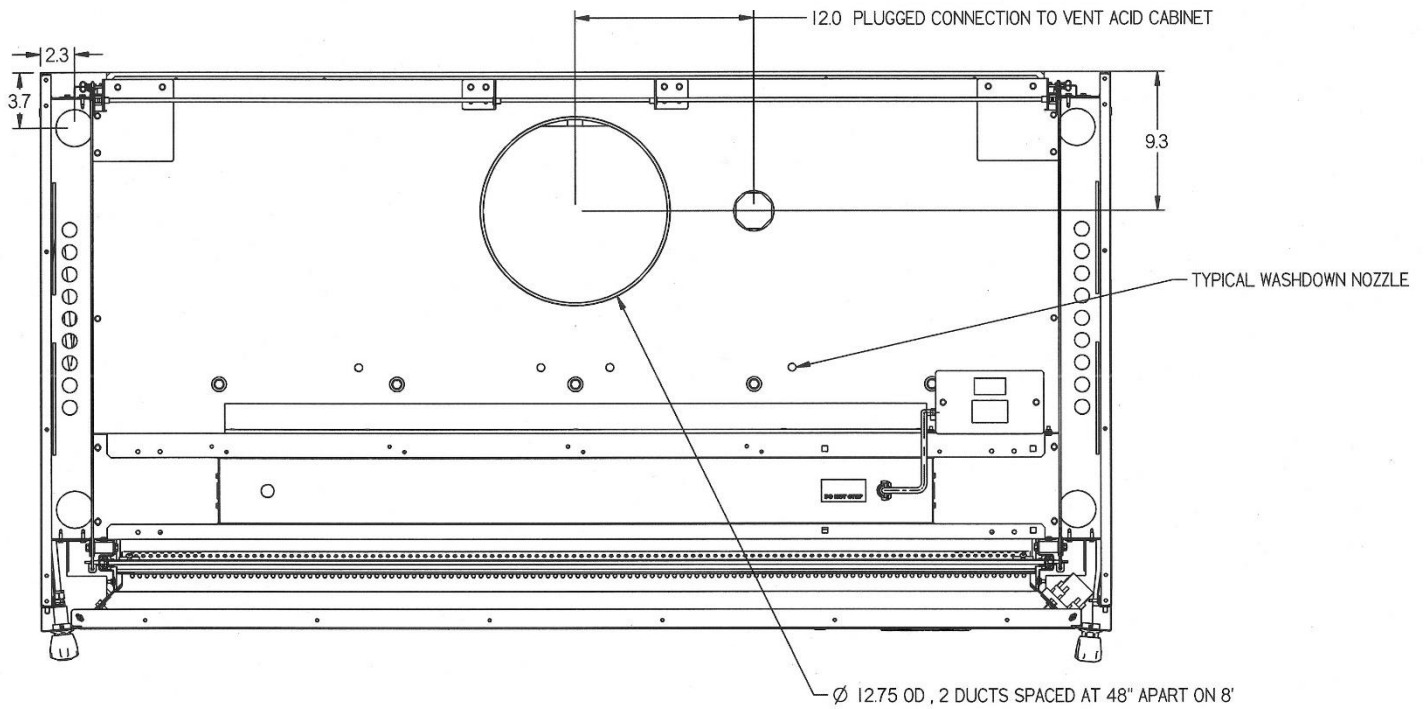
Figure B-3

Appendix B: Protector PVC Dimensions**Figure B-4**

APPENDIX C

PROTECTOR PVC

SPECIFICATIONS

Environmental Conditions


- Indoor use only.
- Maximum altitude: 10,000 feet (3,048 meters).
- Ambient temperature range: 41° to 104°F (5° to 40°C).
- Maximum relative humidity: 80% for temperatures up to 88°F (31°C), decreasing linearly to 50% relative humidity at 104°F (40°C).
- Main supply voltage fluctuations not to exceed $\pm 10\%$ of the nominal voltage.
- Transient over-voltages according to Installation Categories II (Over-voltage Categories per IEC 1010). Temporary voltage spikes on the AC input line that may be as high as 1500V for 115V models and 2500V for 230V models are allowed.
- Used in an environment of Pollution degrees 2 (i.e., where normally only non-conductive atmospheres are present). Occasionally, however, a temporary conductivity caused by condensation must be expected, in accordance with IEC 664.
- Electrical Ratings

<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AMP*</u>
115	1	50/60	10
115	1	50	10
115	1	60	10
230	1	50/60	5
230	1	50	5
230	1	60	5

*Does not include current rating of receptacles.

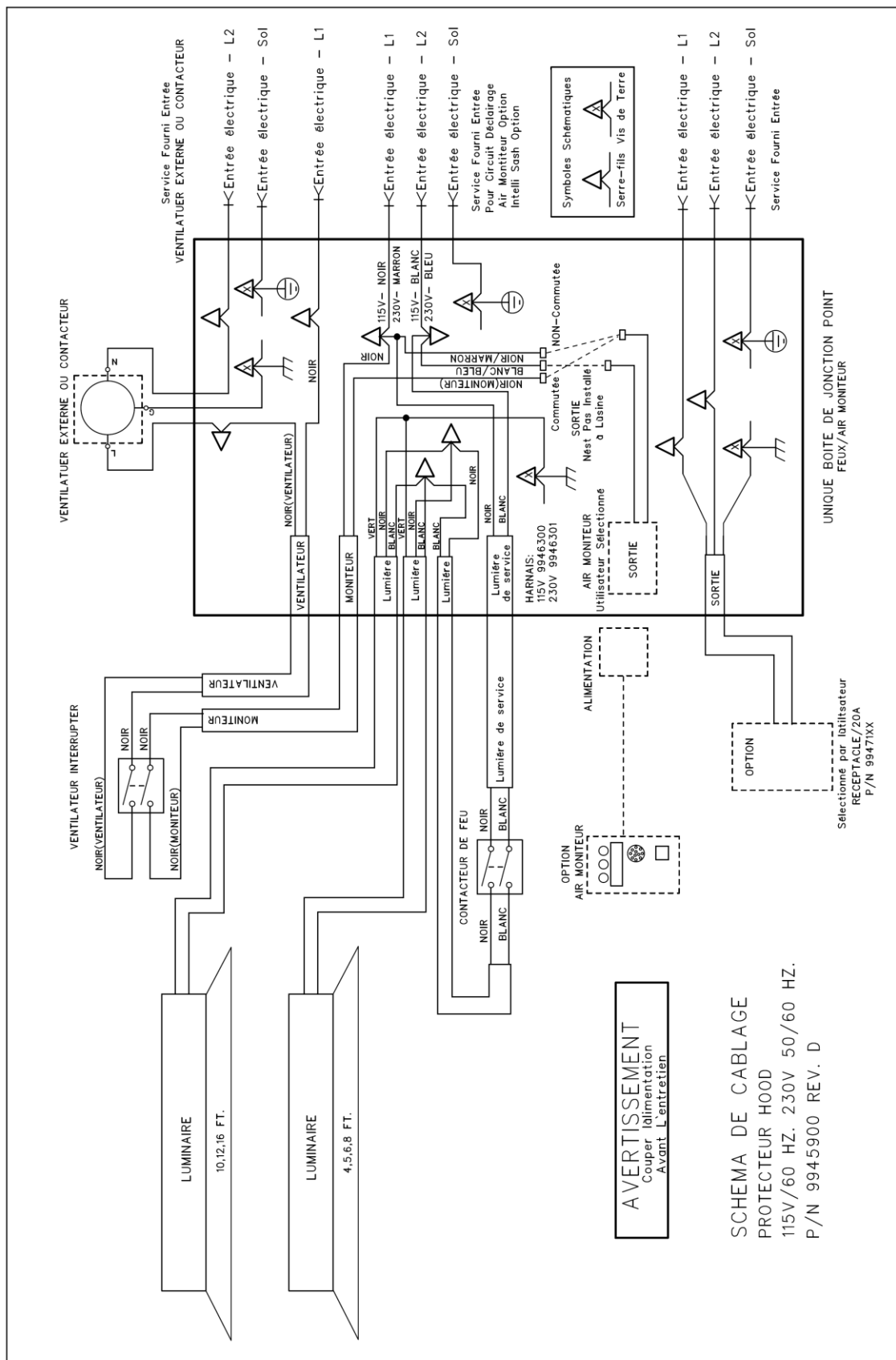Appendix C: Protector PVC Specifications

APPENDIX D

SERIAL NUMBER TAG

DESCRIPTION

Serial tag includes standard information with the following changes:

The portion of the tag for electrical information is labeled with the following units: Volts AC, Phase, Cycle, AMP.

Each hood will include one of the following:

<u>Volts AC</u>	<u>Phase</u>	<u>Cycle</u>	<u>AM</u>  <u>X</u>
115	1	50/60	10X
115	1	50	10X
115	1	60	10X
230	1	50/60	5X
230	1	50	5X
230	1	60	5X

Where  may be a separate label indicating: Caution, See Manual.

The X above will be one of several possible alpha characters as defined in the manual as follows:

- X: No additional duplexes on hood, noted amperage applies.
- B: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- C: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.

- D: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- E: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 115V, single phase, 60Hz, 20 Amps. Each duplex can be wired on a dedicated circuit rated at 20A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 20A electrical service they are wired to.
- F: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- G: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- H: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- J: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **60Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- K: This fume hood includes **one** additional electrical receptacle, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- L: This fume hood includes **two** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

Appendix D: Serial Number Tag Description

- M: This fume hood includes **three** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.
- N: This fume hood includes **four** additional electrical receptacles, individually wired to the field wired box and individually rated as 230V, single phase, **50Hz**, 10 Amps. Each duplex can be wired on a dedicated circuit rated at 10A or the duplexes may be ganged together on the same circuit with the total load of the ganged duplexes not exceeding the 10A electrical service they are wired to.

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APPENDIX E

REFERENCES

Many excellent reference texts and booklets are currently available. The following is a brief listing:

Laboratory Ventilation Standards

Federal Register 29 CFR Part 1910

Non-mandatory recommendations from "Prudent Practices".

- Fume hoods should have a continuous monitoring device
- Face velocities should be between 60-100 linear feet per minute (lfpm)
- Average 2.5 linear feet of hood space per person

Occupational Health and Safety

U.S. Department of Labor

200 Constitution Avenue N.W.

Washington, DC 20210

(202) 523-1452

Industrial Ventilation-ACGIH

- Fume hood face velocities between 60-100 lfpm
- Maximum of 125 lfpm for radioisotope hoods
- Duct velocities of 1000-2000 fpm for vapors, gasses and smoke
- Stack discharge height 1.3-2.0 x building height
- Well designed fume hood containment loss, <0.10 ppm

Industrial Ventilation, A Manual of Recommended Practice.

24th Edition, 2001

American Conference of Governmental Industrial Hygienists

1330 Kemper Meadow drive

Cincinnati, OH 45240-1634

(513) 742-2020

ASHRAE 110-1995 Method of Testing Performance of Fume Hoods

Evaluates fume hood's containment characteristics

- Three part test: Smoke generation, Face velocity profile, Tracer gas release @ 4 liters per minute
- Rated As Manufactured (AM), As Installed (AI) and As Used (AU)

American Society of Heating, Refrigerating, and Air Conditioning Engineers

1791 Tullie Circle N.E.

Atlanta, GA 30329

(404) 636-8400

ANSI Z9.5-2011 Laboratory Standard

Covers entire laboratory ventilation system.

- New and remodeled hoods shall have a monitoring device
- Ductless hoods should only be used with non-hazardous materials

American Industrial Hygiene Association

2700 Prosperity Avenue, Suite 250

Fairfax, VA 22031

(703) 849-8888

SEFA 1-2002

- Fume hood face velocities based on toxicity levels of chemicals
 - Class A – 125 to 150 fpm
 - Class B – 80 to 100 fpm
 - Class C – 75 to 80 fpm
- Test method – face velocity profile and smoke generation

Scientific Equipment & Furniture Association

1028 Duchess Drive

McLean, VA 22102

(703) 538-6007

NFPA 45 – 2002 Fire Protection for Laboratories Using Chemicals

- Laboratory hoods should not be relied on for explosion protection
- Exhaust air from fume hoods should not be recirculated
- Services should be external to the hood
- Canopy hoods only for non-hazardous applications
- Materials of construction should have flame spread of 25 or less
- 80 to 120 fpm to prevent escape

NFPA 30 – 2000 Flammable and Combustible Liquids Code

- Approved cabinets may be metal or wood
- Vent location on cabinets are required
- Venting of cabinets not a requirement

National Fire Protection Association

1 Batterymarch Park

P.O. Box 9101

Quincy, MA 02269-9101

(800) 344-3555

General References

American Conference of Governmental Industrial Hygienists. *Industrial Ventilation, A Manual of Recommended Practice*, Cincinnati, OH

ASHRAE Standard Committee. *ASHRAE Standard* Atlanta: ASHRAE Publications Sales Department, 1995

British Standards Institution, *Laboratory Fume Cupboards*. Parts 1, 2 and 3, London: 1990

Appendix E: References

Department of Labor, Occupational Safety and Health Administration, *29 CFR Part 1910, Occupational Exposures to Hazardous Chemicals in Laboratories, Final Rule*. Vol. 55, No. 21. Washington D.C.:1990

DiBerardinis, L. et al. *Guides for Laboratory Design, Health and Safety Considerations*. Wiley & Sons, 1987

McDermott, Henry, *Handbook of Ventilation for Contaminant Control*, 2nd Edition. Butterworth Publishers, 1985.

Miller, Brinton M. et al. *Laboratory Safety: Principles and Practices*. American Society for Microbiology, Washington, D.C.: 1986

NIH Guidelines for the Laboratory Use of Chemical Carcinogens. NIH Publication No. 81-2385.

Rayburn, Stephen R. *The Foundation of Laboratory Safety, A Guide for the Biomedical Laboratory*. Springer-Verlag, New York: 1990

Sax, N. Irving and Lewis, JR., Richard J. *Rapid Guide to Hazardous Chemicals in the Workplace*. Van Nostrand Reinhold, 1987.

Schilt, Alfred A. *Perchloric Acid and Perchlorates*. The G. Frederick Smith Chemical Company, Columbus, OH: 1979.

Steere, Norman. *CRC Handbook of Laboratory Safety*, 2nd Edition. CRC Press, 1971.

APPENDIX F

VAV OPEN AREA

GUIDELINES

When integrating a Labconco fume hood into a variable air volume (VAV) mechanical system, properly assessing the area opening of the fume hood is critical to ensure correct controller calibration for maintaining face velocity across all sash positions. Depending on your brand of VAV controller, you may be prompted to enter different information to arrive at this correct area.

To simplify this requirement, Labconco offers the following step-by-step procedure to ensure you have the correct information.

1. Locate the serial tag for your fume hood; usually on the lower right of the sash glass.
2. Move to the section of the document below specific to your model number.

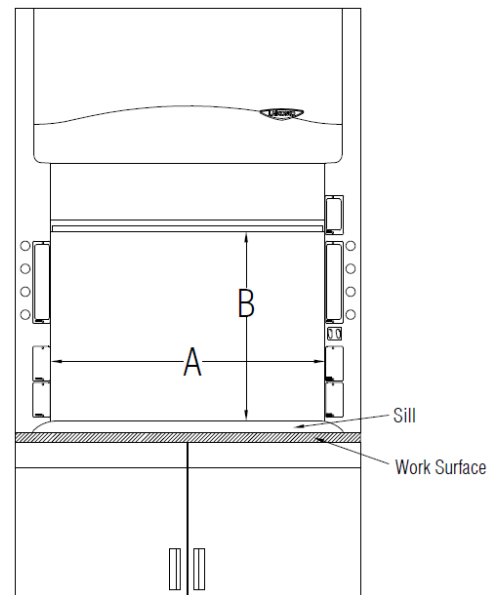
Model numbers beginning 140, 141

1. Determine the free opening area
 - a. Width
 - i. Measure the interior width of the fume hood just behind the plane of the sash glass and enter that here:

Width Opening _____
inches = A

- b. Height*
 - i. Measure the vertical distance between the top of the lower airfoil sill and the bottom of the sash handle, and enter that here:

Vertical Opening _____ inches = B

Appendix F: VAV Open Area Guidelines

c. Free opening area

- i. Multiply the width opening times the vertical opening (A x B), and enter that here:

Free opening area = _____ square inches = C (sq.in)

- ii. Divide this by 144 to arrive at square feet, and enter that here:

Free opening area = _____ square feet = C (sq.ft)

2. Determining bypass or “leakage” area

- a. This is the opening width (A) times 0.34 inches (A x 0.34 = bypass area).

Bypass area = _____ square inches = D (sq in)

- b. Divide this by 144 to arrive at square feet

Bypass area = _____ square feet = D (sq ft)

3. Determining Total Effective Opening Area

- a. This is the Free Opening Area plus the Bypass Opening Area (C + D = Total Effective Opening Area)

Total Effective Area = _____ square inches

Total Effective Area = _____ square feet

* Industry standards call for sash height to be measured from the work surface, however, for the purposes of determining free opening height, measurement is to be taken from the top of the lower airfoil sill to the bottom of the sash handle.